RAYPA

Leading
Lab Technologies

Análisis alimentario
Food analysis
Analyse alimentaire

RAYPA

MÉTODOS DE ANÁLISIS ALIMENTARIO / FOOD ANALYSIS METHODS / MÉTHODES D'ANALYSE ALIMENTAIRE

Kjeldahl

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Kjeldahl

The Kjeldahl method is a chemical analysis process to determine nitrogen content. It is widely used as a reference method to estimate protein content in food.

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Randall/Twisselman

The Randall/Twisselman methods are used to extract substances from a solid sample, usually of lipidic nature, through a solvent.

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Weende

The Weende method is commonly used to quantify crude fibre in a food sample after an acidic and alkaline hydrolysis treatment.

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Van Soest

The Van Soest method is used to determine acid detergent fibre (ADF) and neutral detergent fibre (NDF) content in a sample based on the indigestible residue obtained.

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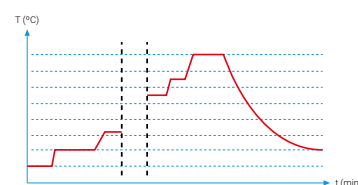
Intended use

- Analysis of proteic nitrogen in organic substances such as nourishing products, fertilizers, seeds, plants, beer, meat, sauces, water or petroleum. See page 3.
- Sample pretreatment (hydrolysis) for fat extraction.
- Soil digestion for metal analysis.

Advantages

- Programmable autostart.
- Built-in temperature regulator which has 9 programs with 18 segments of temperature, ramp and time each.

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- Test tube rack with support (necessary tubes included).
- Fume collector with support.

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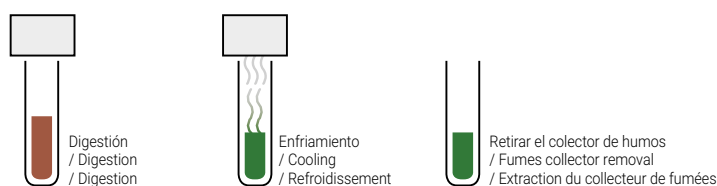
Features

- Quick and safe manipulation of samples.
- Integrated support for sample tubes enabling an easier cooling and saves space.
- Upper chassis made of aluminum alloy and Halar® coating. Metallic bottom chassis painted with epoxy resin. Metal block made of aluminum alloy.
- Thermal insulation: rock wool and ceramic fiber.
- Safety thermostat to prevent overheating.
- Contains calibration port for external probes.
- Ideal for use in combination with the **SCRUBBER Fumes neutralization unit**, (see page 6). Both systems are connected by a flexible tube that joins the Compact digestion system fume collector (MBC) and the Fumes neutralization unit (Scrubber).

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Intended use

- Extraction and neutralization of harmful fumes.

Features

- The gas produced during the digestion is extracted by the action of a vacuum pump. First, the gas undergoes a phase of condensation that acts as preliminary extractor for steams and dragged liquids, avoiding warming or volume increase in posterior wash-up solution.
- The acid or alkaline vapors are washed and neutralized in the next step. In this neutralization phase, most of the unwanted particles are retained through granules of activated charcoal. The used air in the process is conducted to an extractor or to the exterior.
- Case made of stainless steel AISI 304 painted with epoxy resin. Glass objects made of: Borosilicate 3.3. Pipes made of: PVC (Cristalflex®), silicone and Teflon®. Pump made of: PPS, EPDM and Viton®.
- Ideal to use together with **MBC Compact digestion system** (see page 4). Both systems are connected by a flexible tube that joins the Compact digestion system fume collector (MBC) and the Fumes neutralization unit (Scrubber).

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Advantages

- Vacuum pump with acoustic insulation of 48l/min.
- Filtration and neutralization of harmful fumes by a filter of activated carbon.
- Absorption valve manually adjustable from 10 mbar to 800 mbar.
- Adjustable steam suction speed.
- Closed fumes evacuation system.

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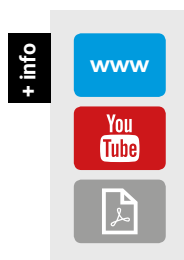
Intended use

• Distillations for the Kjeldahl analysis, ammoniacal nitrogen, alcoholic strength by volume, phenols, volatile acids and cyanides.

Features

- System control by microprocessor with LCD screen.
- Additionally, the system has fixed programs:
 - Ammonium sulphate test.
 - Preheating of the system.
 - Sample waste washing.
- Adaptable system for distillation tubes of different diameters.

Advantages

- Kit for automatic titration included for DNP-3000 and optional for DNP-2000MP.
- Programs: 11 predefined (editable) and 9 free.
- Calibrable pumps by the user.
- Steam generator with water level control.
- Sample suction (only for DNP-2000MP/3000).
- Boric acid pump (only for DNP-2000MP/3000).
- Results in N or total protein (only for DNP-3000).

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Safety

- Open door sensor.
- Distillation tube detection.
- Over-temperature thermostat.
- Overpressure pressure switch.
- Pressure regulator at the cooling water inlet.

Performance

- Measurement range: from 0.1 to 200 mg of N.
- Nitrogen recovery: $\geq 99.5\%$.
- Reproducibility: $\pm 1\%$.
- Distillation time: 4 min/100 ml.
- Electric power: 2000 W (100%).

Programmable parameters and values

- Water for dilution: 0-240 ml.
- NaOH solution: 0-240 ml.
- H_3BO_3 solution: 0-240 ml (only for DNP-2000MP/3000).
- Delay time: 0-30 min.
- Distillation time: 0-30 min.
- Steam flow regulation: 30-100%.

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- 1 distillation tube Ø42 x 300 mm.
- 2 fast connection H_2O and NaOH containers of 10 litres.
- 1 fast connection H_3BO_3 container of 5 litres (only for DNP-2000MP/3000).
- Additional pipe without filter for steam addition (only for DNP-2000MP/3000).
- Automatic titrator (only for DNP-3000).
- Several connection hoses.
- Anti-drip tray.

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Intended use

- Quantification of nitrogen and protein.

It includes

- Magnetic stirring, reaction vessel and complete titrator.

Features

- Potentiometric titrator.
- Multiple titration programs.
- pH calibration.
- Burette autocalibration.
- Results in N or total protein.
- 2 USB ports:
 - 1 to copy tests and extract data.
 - 1 to connect printer, keyboard or mouse.
- Ethernet port: optional software connection.

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Referencia: **SOFT-TITRA**

PC CONNECTION SOFT-TITRA

- Communication software between the titrator and the PC.
- Connection via Ethernet.
- Control to start and stop analysis.
- Show real-time data directly from the workstation.
- Manage data stored locally or on a server (search, compare, delete, print, etc.).
- Export data.

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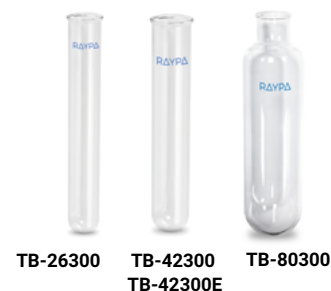
Referencia: **ADAP-BU**

DNP ADAPTER FOR BÜCHI® TUBES

- Adapter to use the Büchi® tubes in the DNP distiller.

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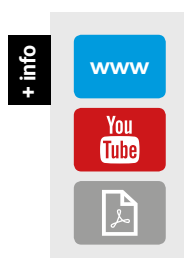
Intended use

• Extraction system for fat or other soluble materials according to Randall or Twisselman methods, based on the evolution of the traditional Soxhlet method. For the analysis of food, feed, soil, detergents, polymers, paper pulp, fibers, gums, textiles, pharmaceuticals and petrochemicals.

Features

- Heating by an electric plate with armoured heater to assure homogeneous distribution of heat.
- Control of temperature by P.I.D. microprocessor action and probe Pt100 class A.
- Over temperature safety control by independent thermocouple.
- Vitón® sealing.
- In accordance with standardized international methods: AOAC, ISO, AACC.
- Case made of stainless steel AISI 304 painted with epoxy resin.

Advantages

- Independent programming of temperature and extraction times for "boiling", "rinsing" and "recovery" phases.

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Technical Data

- Temperature range: room temp. +5 °C to 220 °C.
- Solvent volume: up to 50 ml.
- Solvent recovery: up to 95%.
- Reproducibility: ±1%.
- Extraction thimbles: Øext 26 x 60 mm.
- Programmable boiling extraction time.
- Programmable rinsing extraction time.
- Programmable solvent recovery time.
- Grade of protection: IP65.

Components supplied with the equipment

- 12 extraction vessels + magnetic adapters.
- 6 thimble supports.
- 2 racks for extraction cartridges.
- 2 racks for extraction vessels.
- Aligning rack.
- Insertion handle.
- 2 silicone tubes 1,5m (Ø8x14).
- Cellulose thimbles box (25 units).
- Vessel manipulation rack.
- 6 EPDM seal set.

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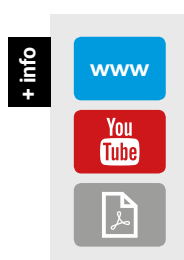
Intended use

• Determination of raw/crude fibre content by *Weende*, acid detergent fibre (ADF) and neutral detergent fibre (NDF) by *Van Soest* and acid detergent lignin (ADL).

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Features

- Extraction and filtration without sample transfer.
- High reproducibility of results.
- Up to 36 analysis/day (*Weende* method).
- Versatile use: samples can be dried/weighted at every phase if needed.
- System for 6 simultaneous samples.
- All electric and mechanic elements are conveniently protected.
- Case made of stainless steel AISI 304 painted with epoxy resin.
- Additionally, it can be used in combination with the **Cold fat extraction system EF-6P** (see page 18), to guarantee the optimum working conditions of the system.

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Technical Data

- Range of measurement: 0.1-100%.
- Reproducibility: $\pm 1\%$ for samples with a fiber content ranging from 1-30%.

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Components supplied with the equipment

- 6 glass crucibles porosity P-2.
- Crucible rack.
- Heater lid.
- Crucible handling support.
- Crucible vessel tong.
- Connection hoses.

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Pressure contribution

- F-6P has a peristaltic pump that provides the necessary conditions for filtering.

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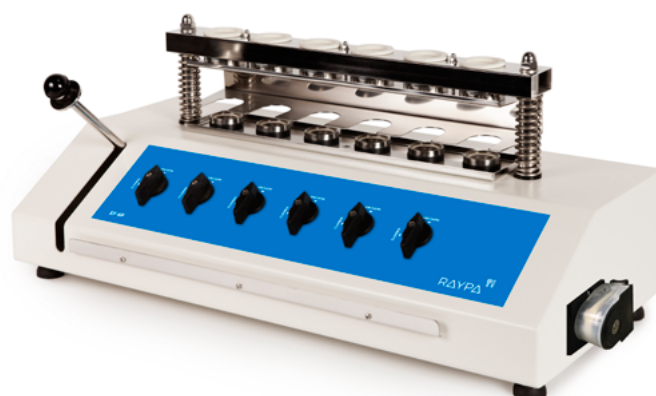
Intended use

- Designed for cold fat extraction with acetone or other solvents.
- Preliminary degrease before hot fibre extraction of samples with more than 1% fat content.

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Features

- Up to 6 simultaneous samples per cycle.
- Case made of stainless steel AISI 304 painted with epoxy resin.
- Ideal for use in combination with the **Manual fibre determination system F-6P** (see page 16).

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Pressure contribution

- EF-6P has a peristaltic pump that provides the necessary conditions for filtering.

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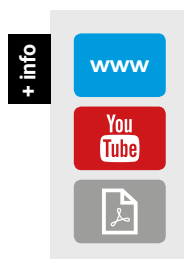
Intended use

• Determination of the alcoholic strength by volume in wines, musts, beers and spirit drinks of high graduation according to the regulation OIV, CEE 2676/90 and CE 2870/2000. Volatile Acidity and Sorbic Acid can also be measured.

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Features

- Distillation by steam dragging.
- Steam generator automatically fed with distilled water, having electronic level control and automatic stop in the case of lack of water in the container.
- Safety thermostat.
- Automatic distillation stop.
- Manual levelling of the distillate avoiding adding distilled water.
- Automatic stop of refrigerating water at the end of distillation.
- Optical display of minimal level of water in the steam generator.
- Button to add distillate manually.
- Case made of stainless steel AISI 304 painted with epoxy resin.

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Technical data

- Distilled water reservoir: 10 l.
- Approx. distillation time: 7 min./200 ml.

Components supplied with the equipment

- 1 distillation tube for maximum sample volume of 100 ml.
- 1 distillation tube for maximum sample volume of 250 ml.
- 1 volumetric flask of 100 ml.
- 1 volumetric flask of 200 ml.
- 1 volumetric flask of 250 ml.
- 1 fast connection container of 10 litres for H₂O.
- Connection hoses.

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TB-42300 TB-80300

GRADILLA 6 TUBOS / 6 TUBES RACK / PORTOIR 6 TUBES

Referencia Reference Référence	Ø x H mm
GRA-640	Ø 42 x 300
GRA-680	Ø 80 x 300

GRA

MATRAZ AFORADO / VOLUMETRIC FLASK / FIOLE JAUGÉE

Referencia / Reference / Référence	Volumen / Volume / Volume ml	Ø x H mm
MA-100	100	Ø 61 x 170
MA-200	200	Ø 75 x 200
MA-250	250	Ø 80 x 220

MA

ALCOHOLÍMETRO / ALCOHOLOMETER / ALCOOLTEST

Referencia Reference Référence	Grado alcohólico Alcoholic strength by volume Titre alcoométrique	Resolución Resolution Résolution
ALC-0-10	0 - 10 %	0,1%
ALC-10-20	10 - 20 %	0,1%
ALC-20-30	20 - 30 %	0,1%
ALC-30-40	30 - 40 %	0,1%
ALC-40-50	40 - 50 %	0,1%
ALC-50-60	50 - 60 %	0,1%
ALC-60-70	60 - 70 %	0,1%
ALC-70-80	70 - 80 %	0,1%
ALC-80-90	80 - 90 %	0,1%
ALC-90-100	90 - 100 %	0,1%

ALC

Avinguda del Vallès, 322
Pol. Ind. "Els Bellots"
08227 Terrassa (Barcelona), Spain

Tel. +34 937 830 720
raypa@raypa.com
www.raypa.com

R. ESPINAR, S.L.