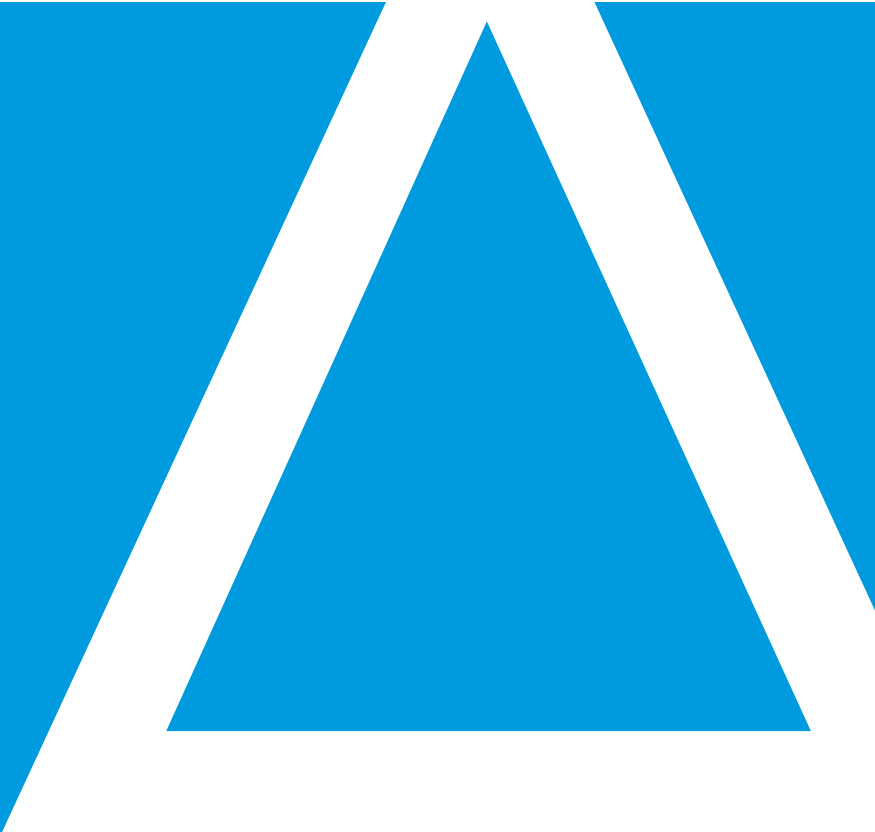
**Leading
Lab Technologies**

CONTACTO
CONTACT
CONTACT

raypa@raypa.com

+34 937 830 720

Avinguda del Vallès, 322
Pol. Ind. "Els Bellots"
08227 Terrassa (Barcelona) Spain

www.raypa.com
www.mediapreparators.com

SOPORTE TÉCNICO
TECHNICAL SERVICE
SERVICE TECHNIQUE

sat@raypa.com
intsat@raypa.com (international)

+34 937 830 720

INDEX

Nuestra empresa	Our company	Notre société	02
Normativas	Regulations	Directives	03

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Sterilization

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04-63

Introducción	Introduction	Introduction	06
Autoclaves verticales	Floor-standing autoclaves	Autoclaves verticaux	15
TOP LINE			
Autoclaves polivalentes	Multipurpose autoclaves	Autoclaves polyvalente	
Autoclaves con sistema de super-secado	Autoclaves with super-drying system	Autoclaves avec système de super-séchage	
Autoclaves con sistema de enfriado rápido	Autoclaves with fast cooling system	Autoclaves avec système de refroidissement rapide	
CLASSIC LINE			
Autoclaves sin secado	Autoclaves without drying	Autoclaves sans séchage	
Autoclaves con secado	Autoclaves with drying	Autoclaves avec séchage	
Autoclaves con prevacíos y secado	Autoclaves with prevacuums and drying	Autoclaves à prévide et séchage	
Autoclaves de sobremesa	Benchtop autoclaves	Autoclaves de paillasse	43
CLASSIC LINE			
Autoclaves sin secado	Autoclaves without drying	Autoclaves sans séchage	
Autoclaves con secado	Autoclaves with drying	Autoclaves avec séchage	
Autoclaves con prevacíos y secado	Autoclaves with prevacuums and drying	Autoclaves à prévide et séchage	
Preparadores de medios	Media preparators	Préparateurs de milieux de culture	58
Incinerador de bacterias	Bacteria incinerator	Incinérateur de bactéries	63

Análisis alimentario

Food analysis

Analyse alimentaire

64-85

Introducción	Introduction	Introduction	66
Sistema de digestión compacto	Compact digestion system	Système compact de digestion	68
Unidad de neutralización de gases	Fumes neutralization unit	Unité de neutralisation de fumées	70
Destiladores nitrógeno-proteína	Protein-nitrogen distiller	Distillateur azote-protéine	72
Sistema de extracción de grasas	Fat extraction system	Système d'extraction de graisse	78
Sistema de determinación manual de fibra	Manual fibre determination system	Système manuel de détermination des fibres	80
Sistema de extracción de grasas en frío	Cold fat extraction system	Système d'extraction de graisse fraîche	82
Destilador enológico	Oenologic distiller	Distillateur oenologique	84

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Incubation and drying

Incubation et séchage

86-103

Introducción	Introduction	Introduction	88
Incubadoras estándar	Standard incubators	Incubateurs standard	92
Incubadoras refrigeradas	Refrigerated incubators	Incubateurs réfrigérées	94
Estufas de secado	Drying ovens	Etuves de séchage	96
Estufas de secado con aire forzado	Forced air drying ovens	Etuves de séchage avec air forcé	98
Estufa de secado al vacío	Vacuum drying oven	Etuve de séchage à vide	100

Baños termostáticos

Thermostatic baths

Bains thermostatiques

104-110

Baños (agua y aceite)	Baths (water & oil)	Bains (eau et huile)	106
Baño de arena	Sand baths	Bains de sable	110

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OUR COMPANY

Raypa, established in 1970, is a company dedicated to the manufacture of laboratory equipment. **Raypa** is specialised in the development, manufacturing, marketing and after-sales service of equipment that requires precise pressure and temperature control. Our main product lines are sterilization equipment, food analysis equipment, incubation and drying equipment and thermostatic baths. All **Raypa** products are manufactured in our Barcelona facilities.

Since its establishment **Raypa** has focused its commercial activity on the international market and currently has a stable distribution network that covers more than 70 countries all over the world.

As a manufacturer that complies with European standards, we have solid procedures in place to ensure the quality and performance of our products. All of them have the **CE** mark for different European quality standards. We are certified by several notified bodies and the company has the UNE-EN-ISO 9001:2015 to guarantee a global quality system.

With more than 40 years of experience, **Raypa** remains faithful to its core values of effort, commitment and service with its customers, suppliers and employees.

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Directiva de aparatos a presión 2014/68/UE.

REGULATIONS

Our instruments comply with the following European directives:

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REGULATIONS

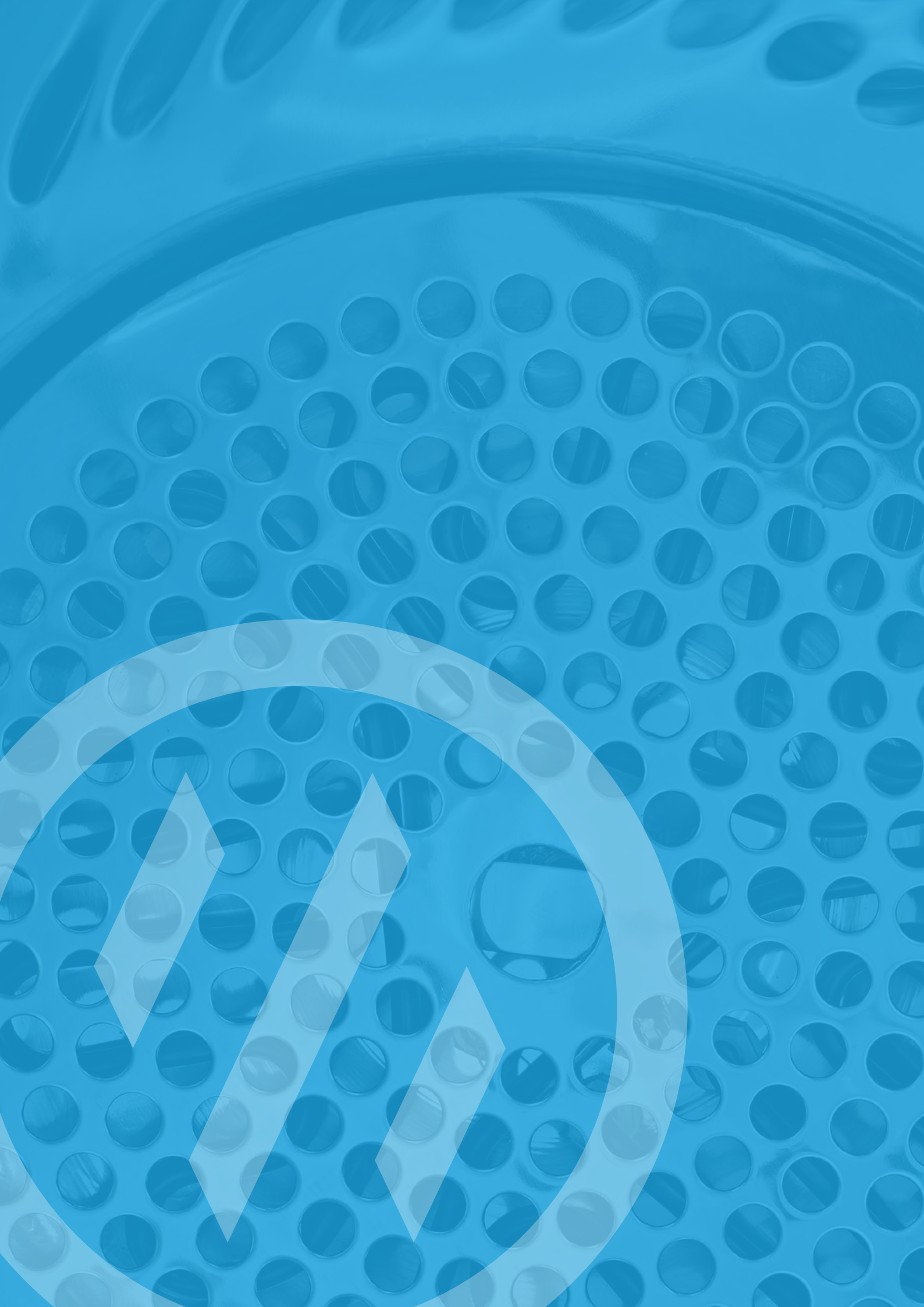
- EN 61010-1 Safety electrical equipment requirements.
- EN 61010-2-010 Particular requirements for heating equipment.
- EN 61010-2-040 Particular requirements for autoclaves using steam.
- EN 61010-2-051 Particular requirements for laboratory equipment for mixing and stirring.
- EN 61326 Electromagnetic compatibility.

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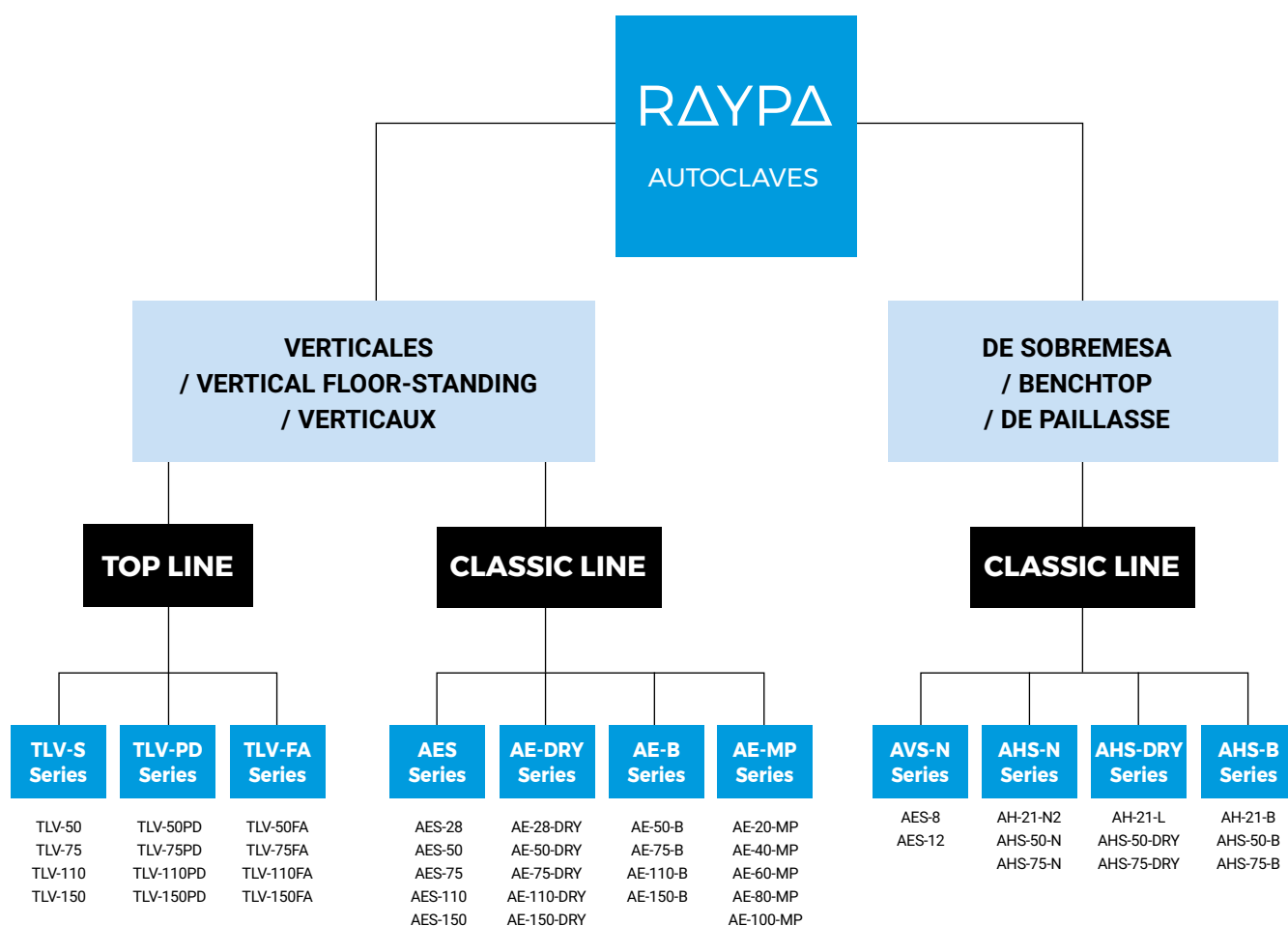
State-of-the-art autoclave technology with **best-in-class connectivity** and multiple process optimization options, designed for research and the biopharmaceutical industry.

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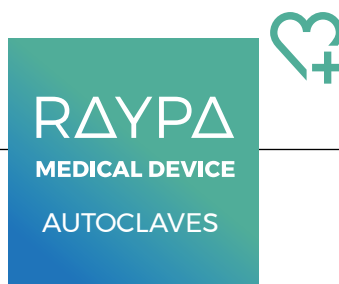
Versatile and **economic** autoclaves for multiple applications, designed for the general laboratory.

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State-of-the-art autoclave technology with **best-in-class connectivity** and multiple process optimization options, designed for biomedical research and the healthcare industry.

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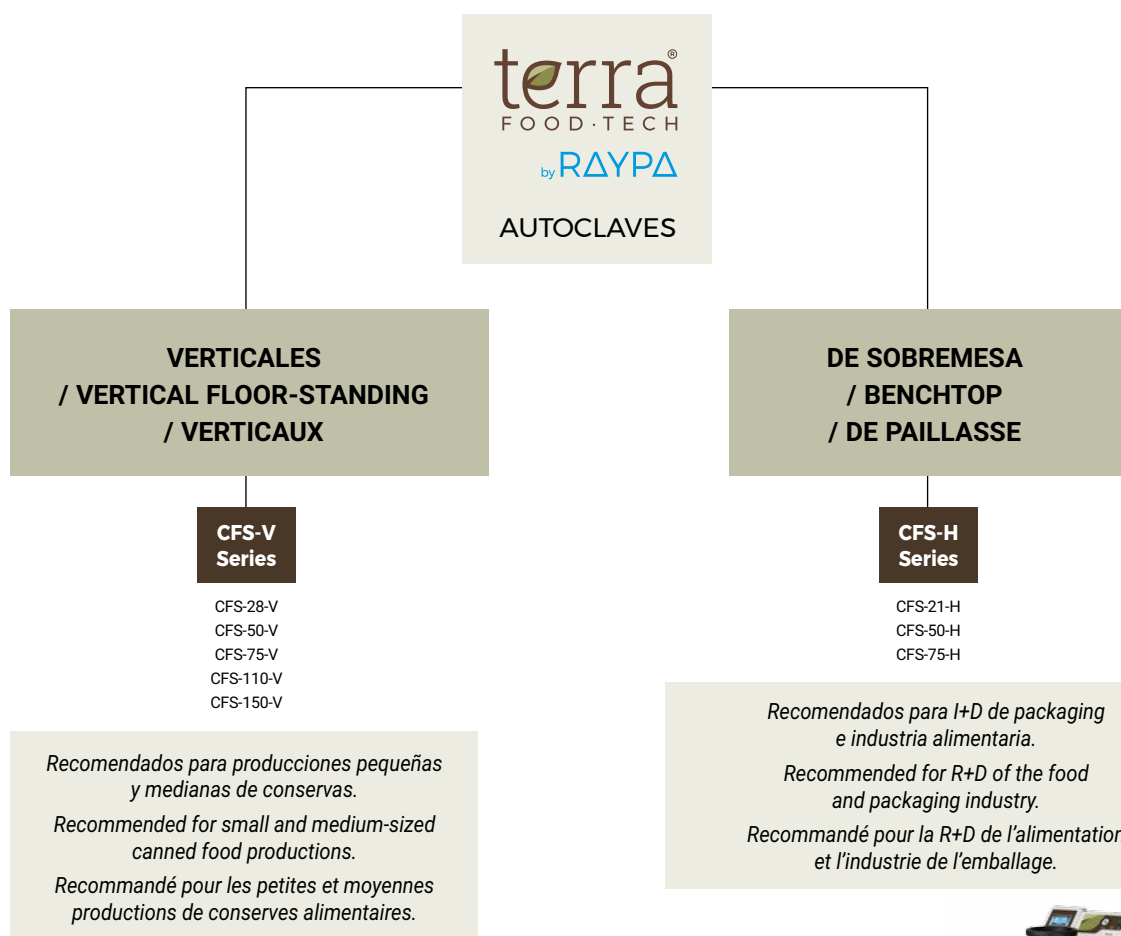
Versatile and **economic** autoclaves for multiple applications, designed for medical laboratories of hospitals and clinics.

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Small retort autoclaves

A new fully integrated autoclave and services brand replacing RAYPA's legacy RFG series, specialised in **small-scale production of canned food** and **food packaging R&D**.

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Main features

- Counterpressure by compressed air system.
- Fast cooling by water shower.
- Sterilization control by ΣF_0 , F_0 , ΣP_0 and P_0
- Extensive in-depth technical support.

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CANNED FOOD AUTOCLAVES

TERRA Food-Tech® allows preserved food sterilization in different types of packaging.

We offer a **wide range of accessories**, customisable according to client needs.

The purchase of an equipment **includes a food consulting service**.

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Type I

AES Series / AVS-N Series / AHS-N Series:

Suitable to sterilize culture medium, glassware, liquids, plastic, metal utensils and laboratory waste bags.

Type II

AE-DRY Series / AHS-DRY Series / TLV-S Series:

Suitable to sterilize unwrapped and wrapped solids, small porous and hollow objects, glassware, plastics, liquids and laboratory waste bags.

Type III

AE-B Series / AHS-B Series / TLV-PD Series:

Indicated for materials that require greater steam penetration to achieve sterilization such as wrapped goods, porous solids, packaged solids, objects with holes and cavities, as well as liquids.

Type IV

AE-MP Series: Indicated for the preparation, sterilization and sterile dispensing of nutrient culture medium for plant tissue culture and microbiology such as liquid agar, peptone water, buffer solutions or lysogenic broth in medium and large scale production settings.

Type V

TLV-FA Series: Multipurpose autoclave with fast cooling system specifically designed for an advanced sterilization of liquids in industrial and research settings of the pharmaceutical and biotechnology industry that sterilizes different types of liquids such as water, solutions and media cultures.

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/ All data on loading capacities of these tables are non-binding guidance to help you choose your autoclave model.

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Differences to consider before choosing a printer

MATRIX SYSTEM

The impression is made by impact, pressing an ink ribbon against the paper. The impression on paper does not fade over time.

THERMAL SYSTEM

Printing is done by transfer of heat on paper. The impression on the paper fades with time.

More information about printers and consumables in the accessories pages.

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TECHNICAL SUMMARY OF OUR VERTICAL FLOOR-STANDING AUTOCLAVES SERIES

		TOP LINE			CLASSIC LINE		
		TLV-S Series	TLV-FA Series	TLV-PD Series	AES Series	AE-DRY Series	AE-B Series
GENERAL CLASSIFICATION	Target client	Research, Biotech, Pharmaceuticals	Research, Biotech, Pharmaceuticals	Research, Biotech, Pharmaceuticals	General laboratory	General laboratory	General laboratory
	Equipment placement	Floor-standing	Floor-standing	Floor-standing	Floor-standing	Floor-standing	Floor-standing
	Load direction	Top-loading	Top-loading	Top-loading	Top-loading	Top-loading	Top-loading
RECOMMENDED TYPE OF LOAD	Liquids and culture media	✓	✓ ✓	✓	✓	✓	✓
	Laboratory waste bags	✓ ✓	✓	✓	✓ ✓	✓	✓
	Porous solids and wrapped loads	-	-	✓ ✓	-	-	✓ ✓
	Biohazardous waste (Red bags)	✓	✓	✓ ✓	-	-	-
	Glassware	✓	✓	✓ ✓	✓	✓ ✓	✓
	Prefilled syringes	-	✓ ✓	-	-	-	-
TRANSFER OF DATA	RS-232	✓	✓	✓	✓	✓	-
	USB & Ethernet	✓	✓	✓	-	-	✓
	Wi-Fi	✓	✓	✓	-	-	-
STERILIZATION CHAMBER AND LID SPECIFICATIONS	Sterilization chamber volume	58 - 166 L	58 - 166 L	58 - 166 L	33 - 175 L	33 - 175 L	55 - 175 L
	External building material	AISI-304	AISI-304	AISI-304	AISI-304	AISI-304	AISI-304
	Sterilization chamber material	AISI-316-L	AISI-316-L	AISI-316-L	AISI-316-L	AISI-316-L	AISI-316-L
	Min. - max. temperature	105 - 140°C	105 - 140°C	105 - 140°C	100 - 134°C	100 - 134°C	105 - 134°C
	Maximum pressure (above atmospheric pressure)	2,6 Barg	2,6 Barg	2,6 Barg	2,1 Barg	2,1 Barg	2,1 Barg
	Automatic locking with pressure	✓	✓	✓	✓	✓	✓
	Mechanism to open the lid	Push-button	Push-button	Push-button	Wheel	Wheel	Wheel
	Direction in which the lid opens	Vertical	Vertical	Vertical	Lateral	Lateral	Lateral
	Mechanically assisted lid opening	✓	✓	✓	-	-	-
	Thermally insulated lid	✓	✓	✓	✓	✓	✓
	Multiple-use water sterilization chamber capacity	-	-	-	2 - 12 L	-	-
WATER MANAGEMENT	Multiple-use water tank capacity	-	-	-	-	9 - 20 L	-
	Single-use water tank capacity	-	-	-	-	-	9 - 20 L
	Automatic water feed from water network	✓	✓	✓	-	0	0
	Water pump to pressurize water network intake	✓	✓	✓	-	0	0
	Water cooled discharge	✓	✓	✓	-	-	-
	Air intake with bacteriological filter	✓	✓	✓	-	✓	✓
OTHER SPECIFICATIONS	Air exhaust with bacteriological filter	✓	✓	✓	-	-	-
	Flexible heart probe	0	✓	0	0	0	0
	Secondary flexible heart probe	0	0	0	-	-	-
	Casters with brakes	✓	✓	✓	0	0	✓
	Pressure gauge	✓	✓	✓	✓	✓	✓
	Electric customisation (115-230M V/230-400T V)	0	0	0	0	0	0
	Special models with augmented power	-	-	-	✓	✓	✓
	Screen display	Touch Panel PC	Touch Panel PC	Touch Panel PC	Digital LCD	Digital LCD	TFT Touchscreen
USER INTERFACE AND MICROPROCESSOR	Screen size	7"	7"	7"	2linex16digits	2linex16digits	5"
	Total number of programs	50	50	50	10	10	50
	Automatic microprocessor control	✓	✓	✓	✓	✓	✓
	Timer start	✓	✓	✓	✓	✓	✓

continued on next page

TECHNICAL SUMMARY OF OUR VERTICAL FLOOR-STANDING AUTOCLAVES SERIES

(continued from)

		TOP LINE			CLASSIC LINE		
		TLV-S Series	TLV-FA Series	TLV-PD Series	AES Series	AE-DRY Series	AE-B Series
STERILIZATION TECHNOLOGY FEATURES	Method to generate steam	Steam generator	Steam generator	Steam generator	Heating elements	Heating elements	Steam generator
	Type of purge	Vacuum	Vacuum	Vacuum	Gravity	Vacuum	Vacuum
	Type of vacuum pump	Membrane	Membrane	Water-ring	-	Membrane	Membrane
	Type of prevacuum	Simple	Simple	Fractioned	-	Simple	Fractioned
	Type of postvacuum	-	-	Fractioned	-	Fractioned	Fractioned
	Heating jacket	-	-	✓	-	✓	✓
	Fast cooling	-	Water coils + Fan	-	-	-	-
	Compressed air system	-	✓	-	-	-	-
	Sterilization controlled by Fo	✓	✓	✓	-	-	-
SPECIAL CYCLES AND PROCESS OPTIMIZATION	Flash cycles (Quick sterilization cycle at high temperatures)	-	-	✓	-	-	✓
	Liquids mode (De-escalated pressurization during cooling phase to avoid liquid boil over)	✓	✓	✓	✓	✓	✓
	Air-over-pressure cycles (Support pressure during cooling phase to avoid liquid loss by evaporation)	-	✓	-	-	-	-
	Agar mode (Temperature holding after cycle ends)	✓	✓	✓	✓	✓	-
	Vacuum drying (Heating jacket + fractioned postvacuum)	-	-	✓	-	✓	✓
	Fast cooling (Faster reduction of temperature)	-	✓	-	-	-	-
	Vacuum leak & Bowie Dick test cycles	-	-	✓	-	-	✓
ADJUSTABLE CYCLE PARAMETERS	Liquids mode	✓	✓	✓	✓	✓	✓
	Agar mode	✓	✓	✓	✓	✓	-
	Temperature of preheating phase	-	-	✓	-	-	-
	Number of prevacuums	✓	✓	✓	-	-	✓
	Temperature and duration of sterilization phase	✓	✓	✓	✓	✓	✓
	Temperature control by heart probe	0	✓	0	0	0	0
	Duration of drying	-	-	✓	-	✓	✓
	Support pressure value during cooling phase	-	✓	-	-	-	-
DATA MANAGEMENT AND ELECTRONIC RECORDS	User accounts and passwords	✓	✓	✓	-	-	-
	Audit trail	0	0	0	-	-	-
	LIMS, Private cloud server	0	0	0	-	-	-
	Internal memory cycle limit	1M	1M	1M	-	-	150 - 200
BATCH PRINTERS	Integrated printer	0	0	0	0	0	0
	External printer	0	0	0	0	0	-
	Label printer and barcode scanner	0	0	0	-	-	-
SERVICES	Third-party qualification (IQ/OQ/PQ)	0	0	0	0	0	0
	Remote live technical assistance	0	0	0	-	-	-
REGULATIONS	ISO 9001	✓	✓	✓	✓	✓	✓
	ISO 13485	0	0	0	0	0	0
	CE marking	✓	✓	✓	✓	✓	✓
	Pressure Equipment Directive 2014/68/EU	✓	✓	✓	✓	✓	✓
	AD 200 Merkblatt Design Codes	✓	✓	✓	✓	✓	✓
	Machine Directive 2006/42/EC	✓	✓	✓	-	-	-
	FDA 21 CFR part 11 & GMP Annex 11 (software and data management conformity)	0	0	0	-	-	-

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Maximize your lab productivity with RAYPAcloud. Thanks to RAYPA, your autoclaves can connect to the internet via Wi-Fi while meeting FDA and GMP requirements for the use of electronically stored data sets and electronic signatures.

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Intended use

• Multipurpose autoclave for solids and liquids designed to cover the sterilization needs in industrial and research settings of the pharmaceutical, chemical and biotechnology industry.

Features

- Sterilization chamber made of AISI-316L stainless steel.
- Max. temperature and pressure: 140°C and 2,6 Barg.
- Integrated steam generator with automatic filling.
- Integrated vacuum pump that allows 1 prevacuum cycle.
- 7" Capacitive color touchscreen.
- FDA & GMP-compliant controller and connectivity to PC, cloud, intranet and LIMS.
- Bacteriological filter for air inlet and outlet.
- Mechanically assisted lid controlled by push-button.
- Automatic water feed from water network.
- Water-cooled discharge.

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Advantages

- Agar mode and programs for liquids.
- Optional main and secondary flexible heart probes to enable sterilization by Fo value.
- Programmable timer start.
- Casters with brakes.
- Unlimited data storage.
- RS-232, USB, Ethernet and Wi-Fi connection.
- Connection to RAYPAcloud.
- Optional integrated or external paper printer, label printer and barcode scanner.
- Compliance with FDA 21 CFR Part 11 and GMP Annex 11.

Safety

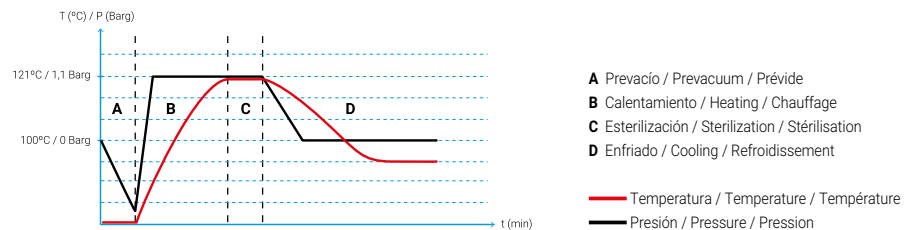
- Safety valve.
- Safety thermostat.
- Lid opening prevention system.
- Open lid sensor.
- Water level detector inside the steam generator.

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Intended use

• Multipurpose autoclave for solids and liquids, with super-drying system for an advanced sterilization of solids in industrial and research settings of the pharmaceutical, chemical and biotechnology industry that sterilize glassware, plastic vials, porous materials and wrapped goods.

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Features

- Sterilization chamber made of AISI-316L stainless steel.
- Maximum temperature and pressure: 140°C and 2,6 Barg.
- Integrated steam generator with automatic filling.
- Super-drying system by heating jacket and liquid-ring pump.
- High performance vacuum pump that allows from 1 to 3 prevacuum cycles.
- 7" Capacitive color touchscreen.
- FDA & GMP-compliant controller and connectivity to PC, cloud, intranet and LIMS.
- Bacteriological filter for air inlet and outlet.
- Automatic water feed from water network.
- Water-cooled discharge.

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Advantages

- Fractioned prevacuum and postvacuum.
- Fast drying function for solids.
- Agar mode and programs for liquids.
- Optional main and secondary flexible heart probes to enable sterilization by Fo value.
- Programmable timer start.
- Casters with brakes.
- Mechanically assisted lid controlled by push-button.
- Unlimited data storage.
- RS-232, USB, Ethernet and Wi-Fi connection.
- Connection to RAYPAcloud.
- Optional integrated or external paper printer, label printer and barcode scanner.
- Compliance with FDA 21 CFR Part 11 and GMP Annex 11.

Safety

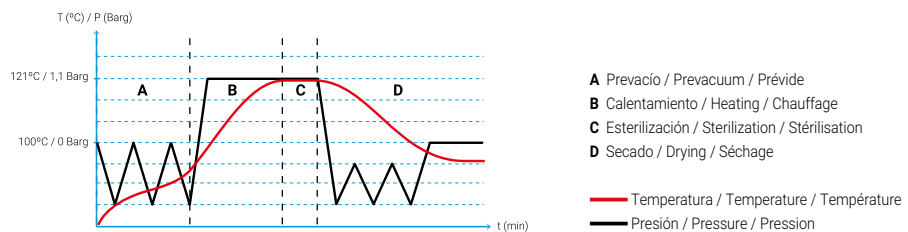
- Safety valve.
- Safety thermostat.
- Lid opening prevention system.
- Open lid sensor.
- Water level detector inside the steam generator.

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Intended use

• Multipurpose autoclave for liquids and solids, with fast cooling system for an advanced sterilization of liquids in industrial and research settings of the pharmaceutical, chemical and biotechnology industry.

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Features

- Sterilization chamber made of AISI-316L stainless steel.
- Max. temperature and pressure: 140°C and 2,6 Barg.
- Integrated steam generator with automatic filling.
- Fast cooling by external water cooling coils and internal radial fan.
- Compressed air system.
- Integrated vacuum pump that allows 1 prevacuum cycle.
- 7" Capacitive color touchscreen.
- FDA & GMP-compliant controller and connectivity to PC, cloud, intranet and LIMS.
- Bacteriological filter for air inlet and outlet.
- Automatic water feed from water network.
- Water-cooled discharge.

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Advantages

- Reduction in the duration of the cooling phase by up to 60% compared to a traditional autoclave.
- Agar mode and programs for liquids.
- Main heart probe included to sterilize by Fo value. Optional secondary heart probe.
- Programmable timer start.
- Casters with brakes.
- Mechanically assisted lid controlled by push-button.
- Unlimited data storage.
- RS-232, USB, Ethernet and Wi-Fi connection.
- Connection to RAYPAcloud.
- Optional integrated or external paper printer, label printer and barcode scanner.
- Compliance with FDA 21 CFR Part 11 and GMP Annex 11.

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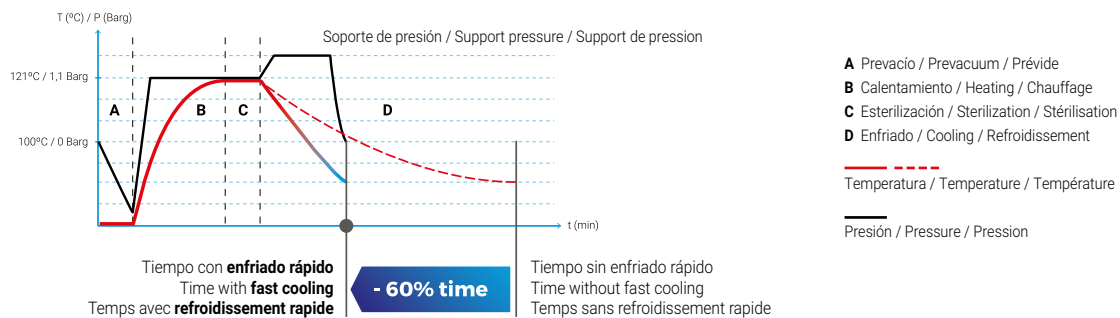
Referencia: **PT-2-TLV**

FLEXIBLE "HEART" PROBE PT-100

- Recommended for sterilizing liquids in flasks of 1 litre or bigger.
- The control is done from the microprocessor.
- Must be installed in our facilities.
- Up to two heart probes can be installed per equipment.

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Consumibles: Papel: **PAPER-ITS**, Tinta: **70934**

DOT MATRIX TABLE TOP PRINTER

- Prints program number, cycle number, temperature, date and hour of the run, and error messages.
- RS-232 connection.

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Consumibles: Papel: **PAPER-IT**, Tinta: **70027**

INTEGRATED THERMAL PRINTER

- Prints program number, cycle number, temperature, date and hour of the run, and error messages.
- Must be installed in our facilities.

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Referencia: **BAR-SCAN**

BARCODE SCANNER

- The barcode scanner allows you to trace the sterilization history of all samples.

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Referencia: **SERVER**

PRIVATE CLOUD SERVER

- Server communicates with the autoclave through the local network of laboratories.
- Automatic backups of all cycle data if RAYPAcloud software is installed.

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Referencia: **FIX-LIFT**

INTEGRATED BASKETS CRANE

- Electrically operated crane made of stainless steel with swivel arm to assist the loading and unloading of heavy items up to 50kg.
- Push-button operation control for ease of use.
- The device is attached to the side of the autoclave.

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Referencia: **MOB-LIFT**

MOBILE BASKETS CRANE

- Electrically operated crane made of stainless steel to assist the loading and unloading of heavy loads up to 50 Kg.
- Push-button operation control for ease of use.
- With swiveling wheels for more maneuverability.

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Referencia: **ECOPUR-500**

ECO-EFFICIENT WATER PURIFIER

• Direct flow eco-efficient water purifier without water accumulation capable of filtering 1.3L/min with LED display user interface control.

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Referencia: **TR-TR**

TRANSPORT TROLLEY

• Auxiliary trolley to assist the loading and unloading of the autoclave.
• Built in chromed iron and plastic.
• The surface of each shelf is textured to prevent the load from shifting.
• Rubber coated wheels to reduce noise.
• Dimensions (LxDxH):
730 x 490 x 700 mm.

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Referencia: **TEST-CT**

STERILIZATION CONTROL TAPE

• Chemical indicator for steam. The change of color indicates that the materials have been processed, without being a guarantee of proper sterilization, Class 1 (to verify that sterilization is effective, additional methods are needed as biological indicators EN ISO 11138).
• Tape roll of 50 m x 19 mm

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Referencia: **TEST-BD**

BOWIE & DICK TEST PACK

• Class B indicator.
• Check the correct steam penetration in the porous loads.
• Indicator printed with non-toxic inks and laminated.
• Box of 20 tests.
For use with the TLV-PD series.

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Intended use

• Suitable to sterilize culture medium, glassware, liquids, plastic, metal utensils and laboratory waste bags.

Features

- Sterilization chamber made of AISI-316L stainless steel. External housing made of AISI-304 stainless steel.
- Equipment controlled by digital PID microprocessor, with 4 predefined and 6 editable programs, adjustable by time, temperature and type of cycle (solids or liquids).
- Alphanumeric LCD screen that shows sterilization parameters, alerts and errors.
- Manual steam release push-button.
- Manual drain valve.
- Manual filling of the sterilization chamber with water.
- Protecting grid for the heater element.
- Cycle time range: 1-250 minutes.

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Advantages

- Agar mode (40 to 80 °C) and programs for liquids.
- Heart probe optional.
- Optional inlet in chamber for an external validation probe.
- Programmable auto-start (24 h).
- Connection to PC.
- RS-232 port to connect to PC or printer.
- Control software optional.
- Optional integrated or external printers.

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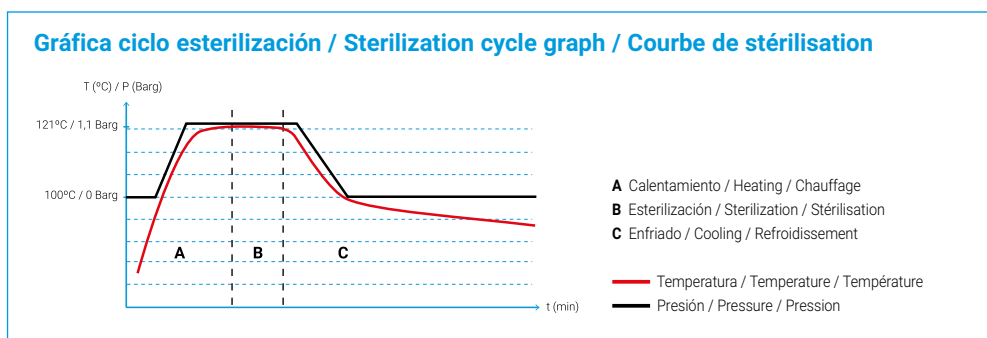
Special models

Available models with increased power to 3200W (AES-28), 6000W (AES-50, AES-75 and AES-110) and 9000W (AES-110 and AES-150). The equipments with increased power reduce the total sterilization cycle time by shortening the heating phase particularly in autoclaves with a high workload.

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Features

- Sterilization chamber made of AISI-316L stainless steel. External housing made of AISI-304 stainless steel.
- Equipment controlled by digital PID microprocessor, with 4 predefined and 6 editable programs, adjustable by time, temperature, drying time and type of cycle (solids or liquids).
- Cycles with initial simple prevacuum and final drying by vacuum pump and heating jacket.
- Alphanumeric LCD screen that shows sterilization parameters, alerts and errors.
- Includes a water tank for an automatic feed of the sterilization chamber. Optional automatic feed of the water tank.
- Protecting grid for the heater element.
- Cycle time range: 1-250 minutes.
- Bacteriological filter for air inlet.

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Safety

- Safety valve and thermostat.
- Hydraulic lid blocking system while positive pressure exists in the chamber.
- Open lid sensor.
- Water level detector in the sterilization chamber.
- Water level detector (min./max.) in the water tank.

Special models

Available models with increased power to 3200W (AE-28DRY), 6000W (AE-50DRY, AE-75DRY and AE-110DRY) and 9000W (AE-110DRY and AE-150DRY). The equipments with increased power reduce the total sterilization cycle time by shortening the heating phase particularly in autoclaves with a high workload.

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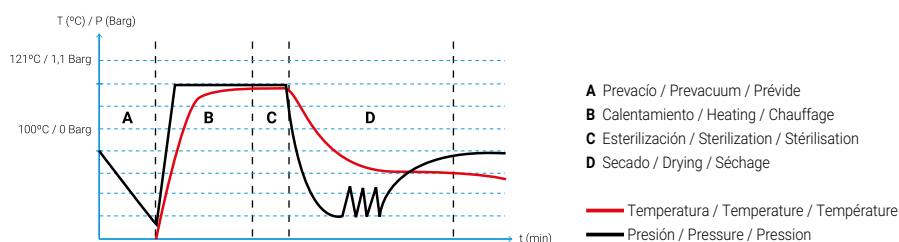
Features

- Sterilization chamber made of AISI-316L stainless steel. External housing made of AISI-304 stainless steel.
- Equipment controlled by digital PID microprocessor, with 12 predefined and 38 editable programs, adjustable by time, temperature, number of prevacuums, drying time and type of cycle (solids or liquids).
- Cycles with fractioned prevacuum (from 1 to 3 prevacuums) and final drying by vacuum pump and heating jacket.
- Immediate production of steam through the built-in steam generator.
- 5" color TFT - LCD touchscreen.
- Includes a single-use water tank for an automatic feed of the sterilization chamber. Optional automatic feed of the water tank.
- Bacteriological filter for air inlet.

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Advantages

- Fractioned prevacuum and automatic drying.
- Programs for liquids.
- Test cycles: Vacuum test and Bowie-Dick test.
- Heart probe (optional).
- Inlet in chamber for an external validation probe.
- Programmable auto-start (date and time).
- Storage of more than 100 cycle data.
- USB port for downloading data and updates. Ethernet port for PC connection.
- Optional integrated printer.

Safety

- Safety valve and thermostat.
- Hydraulic lid blocking system while positive pressure exists in the chamber.
- Open lid sensor.
- Water level detector (min./max.) in the water tank.

Special models

Available models with increased power to 6000W (AE-50B, AE-75B and AE-110B) and 9000W (AE-110B and AE-150B). The equipments with increased power reduce the total sterilization cycle time by shortening the heating phase particularly in autoclaves with a high workload.

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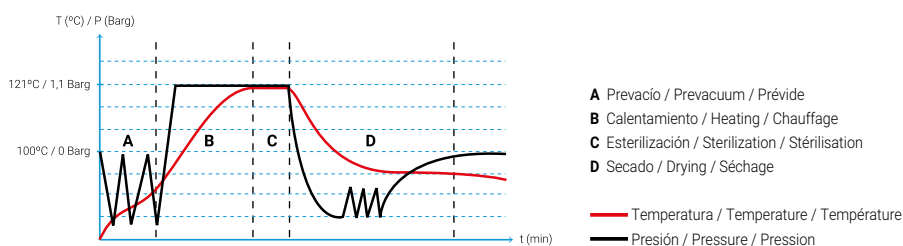
Intended use

- Powerful lift system with adjustable arm to assist the movement of heavy materials into the autoclave.

Features

- Ease of use.
- Compatible with 110L and 150L vertical autoclaves. Consult for other models.
- Up to 200 degrees of rotation.

Safety

- Emergency stop button.
- Motor with auto braking system.

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/ LIQUID COLLECTOR TRAY FOR WIRE BASKETS

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/ UNPERFORATED STAINLESS STEEL BASKETS FOR WASTE STERILIZATION

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/ STAINLESS STEEL "SCHIMMELBUSCH" DRUM FOR MEDICAL INSTRUMENTS

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FLEXIBLE "HEART" PROBE PT-100

- Recommended for sterilizing liquids in flasks of 1 litre or bigger.
- The control is done from the microprocessor.
- Must be installed in our facilities.
- Compatible with AES, AE-DRY and AE-B series.

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Referencia: **EXT-TP**

EXTERNAL PROBE ADAPTER

- External adapter for continuous validation processes that allows the access of an external probe (Ø 3-6 mm) to obtain a temperature reading independent from that of the equipment microprocessor.
- It is located on the lid of the autoclave.

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AUTOMATIC WATER FILLING KIT

- Water pump to automate the feed of the water tank with purified water in AE-DRY series and AE-B series autoclaves.
- Compatible with installations that either have a purified water network, a purified water tank or facilities that have an unpurified water network, in the latter case, the kit should be supplied with two other accessories: water purifier (ECOPUR-500) and purified water tank (TANK-KLL).
- Must be installed in our facilities.

References: **KLL** (AE-DRY series) and **KLL-B** (AE-B series).

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Referencia: **ECOPUR-500**

ECO-EFFICIENT WATER PURIFIER

- Direct flow eco-efficient water purifier without water accumulation capable of filtering 1.3L/min with LED display user interface control.
- The installation of this accessory in the AE-DRY, AE-B and AE-MP autoclaves requires the joint installation of the external tank (TANK-KLL) and the automatic filling system (KLL) corresponding to each model.

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Referencia: **4WHBR**

SET OF 4 WHEELS (2 WITH BRAKES)

- Enables an easy movement of the equipment between different locations. Includes brakes on two of the casters.
- Compatible with AES and AE-DRY series. Already included in AE-B.

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Referencia: **TR-TR**

TRANSPORT TROLLEY

- Auxiliary trolley to assist the loading and unloading of the autoclave.
- Built in chromed iron and plastic.
- The surface of each shelf is textured to prevent the load from shifting.
- Rubber coated wheels to reduce noise.
- Dimensions (LxDxH): 730 x 490 x 700 mm.

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Referencia: **TANK-AE**

CONDENSATES BOTTLE

- 12L bottle with tap to collect condensates during purge phase and also to collect waste water while cleaning the sterilization chamber. Useful in situations where no easy drainage is available.
- For use with the AES series.

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Referencia: **SW8000**

PC CONNECTION SW8000

- Communication software between the equipment and the PC.
- Used with Ethernet connection.
- USB adapter.
- Compatible with AE-B series.

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Referencia: **SW7000**

PC CONNECTION SW7000

- Communication software between the equipment and the PC.
- Used with RS-232 connection.
- USB adapter.
- Compatible with AES and AE-DRY series.

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Referencia: **TEST-CT**

STERILIZATION CONTROL TAPE

- Chemical indicator for steam. The change of color indicates that the materials have been processed, without being a guarantee of proper sterilization, Class 1 (to verify that sterilization is effective, additional methods are needed as biological indicators EN ISO 11138).
- Tape roll of 50 m x 19 mm

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Referencia: **TEST-BD**

BOWIE & DICK TEST PACK

- Class B indicator.
- Check the correct steam penetration in the porous loads.
- Indicator printed with non-toxic inks and laminated.
- Box of 20 tests.
- For use with the AE-B series.

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TECHNICAL SUMMARY OF OUR BENCHTOP AUTOCLAVES SERIES

		CLASSIC LINE							
		AVS-N Series		AHS-N Series		AHS-DRY Series		AHS-B Series	
AVAILABLE MODELS		AES-8	AES-12	AH-21-N2	AHS-50-N AHS-75-N	AH-21-L	AHS-50-DRY AHS-75-DRY	AH-21-B	AHS-50-B AHS-75-B
GENERAL CLASSIFICATION	Target client	General laboratory		General laboratory		General laboratory		General laboratory	
	Equipment placement	Benchtop		Benchtop		Benchtop		Benchtop	
	Load direction	Top-loading		Front-loading		Front-loading		Front-loading	
RECOMMENDED TYPE OF LOAD	Liquids and culture media	-	✓	✓		✓		✓	
	Laboratory waste bags	✓	✓	✓	✓	✓		✓	
	Porous solids and wrapped loads	-		-		-		✓	✓
	Glassware	✓		✓		✓	✓	✓	
TRANSFER OF DATA	RS-232	✓		✓		✓		-	
	USB & Ethernet	-		-		-		✓	
STERILIZATION CHAMBER AND LID SPECIFICATIONS	Sterilization chamber volume	8 L	15 L	22 L	55 - 79 L	22 L	55-79 L	22 L	55 - 79 L
	Sterilization chamber material	AISI-18/10	AISI-316-L	AISI-316-L		AISI-316-L		AISI-316-L	
	Min. - max. temperature	100 - 127°C	100 - 134°C	100 - 134°C		100 - 134°C		105 - 134°C	
	Maximum pressure (Above atmospheric pressure)	1,5 Barg	2,1 Barg	2,1 Barg		2,1 Barg		2,1 Barg	
	Automatic locking with pressure	✓		✓		✓		✓	
	Mechanism to open the lid	Bayonet	Wheel	Handle	Wheel	Handle	Wheel	Handle	Wheel
	Direction in which the lid opens	Vertical		Frontal		Frontal		Frontal	
	Thermally insulated lid	✓		✓		✓		✓	
WATER MANAGEMENT	Multiple-use water sterilization chamber capacity	1,2 L	2 L	-		-		-	
	Multiple-use water tank capacity	-		6 L	10 L	6 L	10 L	2,6 L	6 L
	Single-use water tank capacity	-		-		-		6 L	10 L
	Automatic water feed from water network	-		-		0		0	
	Water pump to pressurize water network intake	-		-		0		0	
OTHER SPECIFICATIONS	Air intake with bacteriological filter	-		-		✓		✓	
	Flexible heart probe	-	0	-	0	0		0	
	Pressure gauge	✓		✓		✓		✓	
	Internal memory cycle data storage limit	-		-		-		150-200	
	Custom electrical features (115-230M V / 230-400T V)	0		0		0		0	
USER INTERFACE AND MICROPROCESSOR	Screen display	Digital LCD		Digital LCD		Digital LCD		TFT Touchscreen	
	Screen size	1line x 3digits	2line x 16digits	1line x 3digits	2line x 16digits	2line x 16digits		5"	
	Total number of programs	1	10	1	10	10		50	
	Automatic microprocessor control	✓		✓		✓		✓	
	Timer start	✓		✓		✓		✓	

continued on next page

TECHNICAL SUMMARY OF OUR BENCHTOP AUTOCLAVES SERIES

(continued from)

		CLASSIC LINE						
		AVS-N Series		AHS-N Series		AHS-DRY Series		AHS-B Series
AVAILABLE MODELS		AES-8	AES-12	AH-21-N2	AHS-50-N AHS-75-N	AH-21-L	AHS-50-DRY AHS-75-DRY	AH-21-B AHS-50-B AHS-75-B
STERILIZATION TECHNOLOGY FEATURES	Method to generate steam	Heating elements		Heating elements		Heating elements		Steam generator
	Type of purge	Gravity		Gravity		Vacuum		Vacuum
	Type of vacuum pump	-		-		Membrane		Membrane
	Type of prevacuum	-		-		Simple		Fractioned
	Type of postvacuum	-		-		Fractioned		Fractioned
	Heating jacket around sterilization chamber	-		-		✓		✓
SPECIAL CYCLES AND PROCESS OPTIMIZATION	Flash cycles (Quick sterilization cycle at high temperatures)	-		-		-		✓
	Liquids mode (De-escalated pressurization during cooling phase to avoid liquid boil over)	✓		✓		✓		✓
	Agar mode (Temperature holding after cycle ends)	-	✓	-	✓	✓		-
	Vacuum drying (Heating jacket + fractioned postvacuum)	-		-		✓		✓
	Vacuum leak & Bowie Dick test cycles	-		-		-		✓
ADJUSTABLE CYCLE PARAMETERS	Liquids mode	-	✓	✓		✓		✓
	Agar mode	-	✓	-	✓	✓		-
	Number of prevacuums	-		-		-		✓
	Temperature and duration of sterilization phase	✓		✓		✓		✓
	Temperature control by heart probe	-	0	-	0	0		0
	Duration of drying	-		-		✓		✓
DATA MANAGEMENT	Internal memory cycle limit	-		-		-		150-200
BATCH PRINTERS	Integrated printer	-	0	-	0	0		0
	External printer	-	0	-	0	0		-
SERVICES	Third-party qualification (IQ/OQ/PQ)	0		0		0		0
REGULATIONS	ISO 9001	-		-		-		-
	ISO 13485	-	0	0		0		0
	CE marking	✓		✓		✓		✓
	Pressure Equipment Directive 2014/68/EU	✓		✓		✓		✓
	AD 200 Merkblatt Design Codes	✓		✓		✓		✓

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Intended use

- Suitable to sterilize culture medium, glassware, liquids, plastic, metal utensils and laboratory waste.

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Features

- Sterilization chamber made of 18/10 stainless steel. External housing made of AISI-304 stainless steel.
- Equipment controlled by digital PID microprocessor, adjustable by time and temperature.
- Alphanumeric LCD screen that shows sterilization parameters.
- RS-232 port to connect to PC.
- Manual valves for drainage and unsteaming.
- Cycle time range: 0 - 500 h.
- Manual filling of the sterilization chamber.
- Protecting grid for the heater element.

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Safety

- Safety valve and thermostat.
- Open lid blocking system.

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Features

- Sterilization chamber made of AISI-316L stainless steel. External housing made of AISI-304 stainless steel.
- Equipment controlled by digital PID microprocessor, with 4 predefined and 6 editable programs, adjustable by time and temperature.
- Alphanumeric LCD screen that shows sterilization parameters, alerts and errors.
- RS-232 port to connect PC or printer.
- Manual filling of the sterilization chamber with water.
- Protecting grid for the heater element.
- Manual steam release push-button.
- Manual valve to drain.
- Cycle time range: 1-250 minutes.

Advantages

- Agar mode (40 to 80 °C) and programs for liquids.
- Optional heart probe.
- Inlet in chamber for an external validation probe.
- Programmable auto-start (24 h).
- Connection to PC.
- Optional control software.
- Optional integrated or external printers.

Safety

- Safety valve and thermostat.
- Hydraulic lid blocking system while positive pressure exists in the chamber.
- Open lid sensor.

Components supplied

- Stainless steel wire basket.
- For AES-8: CV-8 and for AES-12: CV-12

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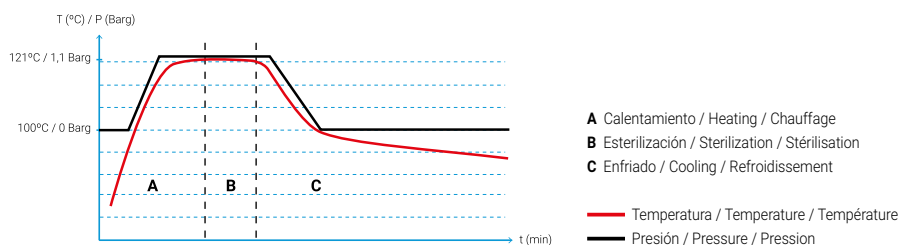
CV
Cesto incluido
/ Basket included
/ Panier inclus

Gráfica ciclo esterilización / Sterilization cycle graph / Courbe de stérilisation

Referencia Reference Référence	Volumen total / útil Total volume / usable Volume total / utile l	Dimensiones útiles cámara Usable chamber dimensions Dimensions utile chambre Ø x H mm	EXTERIOR L x D x H mm	Temp. °C	Presión máx. Max. pressure Pression max. Barg	Potencia Power Puissance W	Peso bruto Gross weight Poids brut Kg	Tensión* Voltage Tension V	Frecuencia Frequency Fréquence Hz
AES-8	8/6,8	220 x 180	410 x 355 x 430	100 - 127	1,5	1000	11	230	50/60
AES-12	15/13	250 x 280	490 x 475 x 630	100 - 134	2,1	1000	35	230	50/60

* Consultar por otras tensiones / Ask for other voltages / Demander d'autres tensions

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Features

- Sterilization chamber made of AISI-316L stainless steel.
- Equipment controlled by digital PID microprocessor, cycles adjustable by time and temperature.
- Alphanumeric LCD screen that shows sterilization parameters.
- RS-232 port for PC.
- Includes a manually-filled water tank that automatically feeds the sterilization chamber.
- Time range: 0-500 h

Advantages

- Switch for solids or liquids cycles.
- Connection to PC.
- Optional control software.
- Programmable auto-start: 500h.

Safety

- Safety valve and thermostat.
- Hydraulic lid blocking system while positive pressure exists in the chamber.
- Open lid sensor.

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Features

- Sterilization chamber made of AISI-316L stainless steel.
- Equipment controlled by digital PID microprocessor with 4 predefined and 6 editable programs, adjustable by time and temperature.
- Alphanumeric LCD screen that shows sterilization parameters, alerts and errors.
- RS-232 port to connect PC or printer.
- Includes a manually-filled water tank that automatically feeds the sterilization chamber.
- Manual steam release push-button.
- Manual drain valve.
- Cycle time range: 1-250 min.

Advantages

- Agar mode (40 to 80 °C) and programs for liquids.
- Optional heart probe.
- Connection to PC.
- Optional control software.
- Programmable auto-start: 24 h
- Optional integrated or external printers.

Safety

- Safety valve and thermostat
- Hydraulic lid blocking system while positive pressure exists in the chamber.
- Open lid sensor.

Components supplied**AH-21-N2**

- Support for 4 trays.
- 3 stainless steel perforated trays.
- Auxiliary tray.
- A holding clamp.
- 1.5 m flexible tube for drainage.

AHS-50-N & AHS-75-N

- Support for 5 trays.
- 2 stainless steel perforated trays.
- Auxiliary tray.
- 1.5 m flexible tube for drainage.

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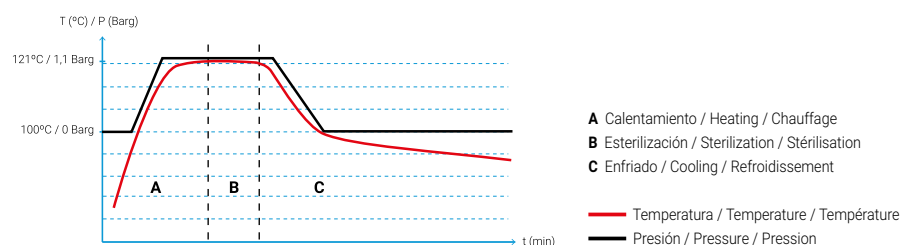
Intended use

• Suitable to sterilize unwrapped and wrapped solids, small porous and hollow objects, glassware, plastics, liquids and laboratory waste bags.

Features

- Sterilization chamber made of AISI-316L stainless steel.
- Equipment controlled by digital PID microprocessor, with 4 predefined and 6 editable programs, adjustable by time, temperature, drying time and type of cycle (solids or liquids).
- Cycles with initial simple prevacuum and final drying by vacuum pump and heating jacket.
- Alphanumeric LCD screen that shows sterilization parameters, alerts and errors.
- RS-232 port to connect PC or printer.
- Includes a water tank that automatically feeds the sterilization chamber. Optional automatic feed of the water tank.
- Cycle time range: 1-250 min.
- Bacteriological filter for air inlet.

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Safety

- Safety valve and thermostat.
- Hydraulic lid blocking system while positive pressure exists in the chamber.
- Open lid sensor.
- Water level detector.

Components supplied**AH-21-L**

- Support for 4 trays.
- 3 stainless steel perforated trays.
- Auxiliary tray.
- Holding clamp.
- 1.5 m flexible tube for drainage.
- Bag support.

AHS-50-DRY & AHS-75-DRY

- Support for 5 trays.
- 2 stainless steel perforated trays.
- Auxiliary tray.
- 1.5 m flexible tube for drainage.

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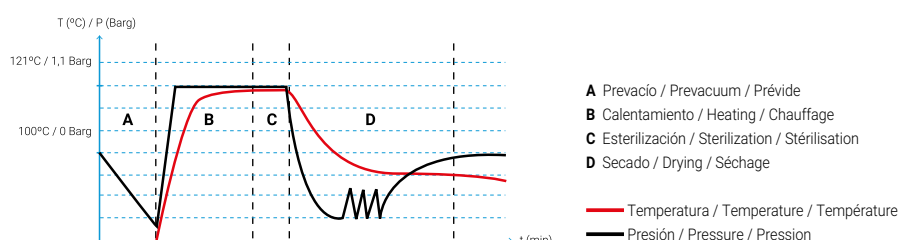
Intended use

• Indicated for materials that require greater steam penetration to achieve sterilization such as wrapped goods, porous solids, packaged solids, objects with holes and cavities, as well as liquids.

Features

- Sterilization chamber made of AISI-316L stainless steel.
- Equipment controlled by digital PID microprocessor, with 12 predefined and 38 editable programs, adjustable by time, temperature, number of prevacuums, drying time and type of cycle (solids or liquids).
- Cycles with fractioned vacuum (1-3 prevacuums) and final drying by vacuum pump and heating jacket.
- Steam production by steam generator.
- Color alphanumeric touchscreen that shows sterilization parameters, alerts and errors.
- Includes clean and used water tanks for an automatic feed of the sterilization chamber. Optional automatic feed of the clean water tank.
- Bacteriological filter for air inlet.

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Advantages

- Fractioned prevacuum and automatic drying.
- Programs for liquids.
- Test cycles: Vacuum Test and Bowie-Dick Test.
- Optional heart probe.
- Connection to PC.
- USB port for downloading data and updates. Ethernet port for PC connection.
- Internal memory for more than 100 cycle data.
- Optional control software.
- Programmable auto-start (date and time).
- Optional integrated printer.

Safety

- Safety valve, thermostat and pressure switch.
- Hydraulic lid blocking system while positive pressure exists in the chamber.
- Open lid sensor.
- Water level detector.

Components supplied**AH-21-B**

- Support for 4 trays.
- 3 stainless steel perforated trays.
- Auxiliary tray.
- Holding clamp.
- 1.5 m flexible tube for drainage.
- Bag support.

AHS-50-B & AHS-75-B

- Support for 5 trays.
- 2 stainless steel perforated trays.
- Auxiliary tray.
- 1.5 m flexible tube for drainage.

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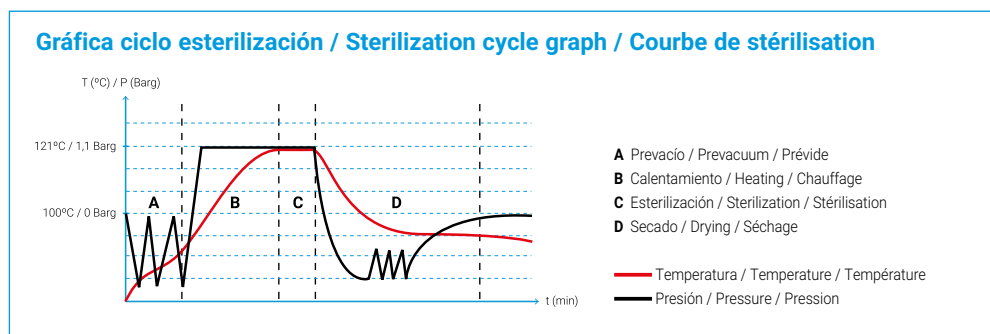
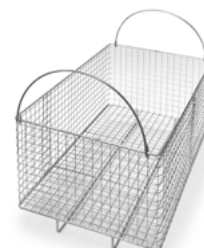
Referencias: **PT-2-AH****FLEXIBLE "HEART"
PROBE PT-100**

- Recommended for sterilizing liquids in flasks of 1 litre or bigger.
- The control is done from the microprocessor.
- Must be installed in our facilities.
- Compatible with AVS-N (AES-12 model), AHS-N (AHS-50-N and AHS-75-N models), AHS-DRY and AE-B series.

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Referencias: **KLL-21 y KLL-AHS****AUTOMATIC WATER
FILLING KIT**

- Water pump to automate the feed of the water tank with purified water in benchtop autoclaves.
- Compatible with installations that either have a purified water network, a purified water tank or facilities that have an unpurified water network, in the latter case, the kit should be supplied with two other accessories: water purifier (ECOPUR-500) and purified water tank (TANK-KLL).
- Must be installed in our facilities.

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Referencia: **DEM-4****WATER DISTILLER**

- Forced air water distiller with stainless steel interior, and a capacity of 4L and a distillation volume output of 1.5L/h.

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Referencia: **SW8000**

PC CONNECTION SW8000

- Communication software between the equipment and the PC.
- Used with Ethernet connection.
- USB adapter.
- Compatible with AHS-B series.

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Referencia: **SW7000**

PC CONNECTION SW7000

- Communication software between the equipment and the PC.
- Used with RS-232 connection.
- USB adapter.
- Compatible with AVS-N, AHS-N and AE-DRY series.

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Referencia: **TEST-CT**

STERILIZATION CONTROL TAPE

- Chemical indicator for steam. The change of color indicates that the materials have been processed, without being a guarantee of proper sterilization, Class 1 (*to verify that sterilization is effective, additional methods are needed as biological indicators EN ISO 11138*).
- Tape roll of 50 m x 19 mm

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Referencia: **TEST-BD**

BOWIE & DICK TEST PACK

- Class B indicator.
- Check the correct steam penetration in the porous loads.
- Indicator printed with non-toxic inks and laminated.
- Box of 20 tests.
- For use with the AHS-B series.

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Intended use

• Preparation, sterilization and sterile dispensing of nutrient culture medium for plant tissue culture and microbiology such as liquid agar, peptone water, buffer solutions or lysogenic broth in medium and large scale production settings.

Features

- Sterilization chamber made of AISI-316L stainless steel.
- Fast cooling by water cooling coils.
- Optional compressed air system.
- 5" Color TFT-LCD touchscreen.
- 50 programs available, adjustable by time, temperature or Fo, with the option of temperature control by flexible heart probe or direct chamber temperature. You can also keep constant the temperature at the end of cycle, adjust the dispensing temperature, the stirring speed and the counterpressure value while dispensing.
- Precise dispensing controlled by pedal or by peristaltic pump pulses for repetitive filling operations. Optional external dispensing control with the DW-MP accessory (page 60).
- Direct feed from water network or optional manual addition in the second sterilization chamber.
- Bacteriological filter for air inlet.

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Advantages

- Integration in a single equipment of the preparation, sterilization, cooling and dispensing.
- Reduction in the time of the cooling phase by up to 60% compared to a traditional autoclave.
- Incorporated heart probe, sterilization controlled by Fo.
- Optional control and registry of pH, with real-time monitoring on screen.
- Powerful magnetic stirrer adjustable from 50 to 200 rpm.
- Programmable peristaltic pump.
- Easy extraction and cleaning of the media chamber thanks to integrated handle grips.
- Dispensing lines can be cleaned each time with steam pulses.
- USB port (for downloading data and charging new updates) and Ethernet port for PC connection.
- Optional integrated printer.

Safety

- Safety valve and thermostat.
- Hydraulic lid blocking system while existing positive pressure in the chamber.
- Open lid sensor.
- Water level detector.

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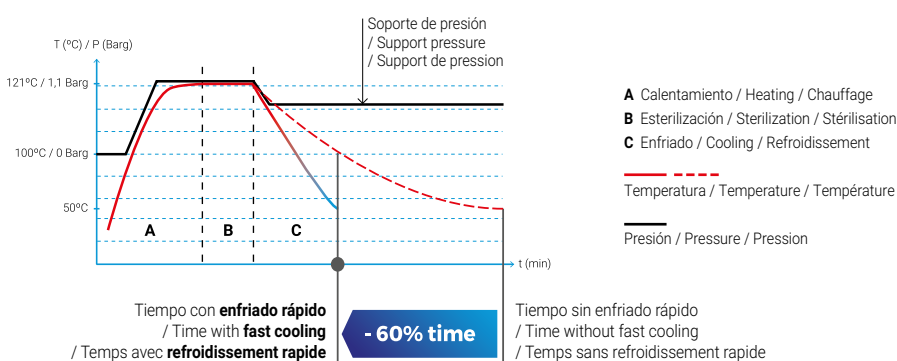
PT-2
Sonda corazón incluida
/ Included heart probe
/ Sonde flexible inclus

Gráfica ciclo esterilización / Sterilization cycle graph / Courbe de stérilisation

Referencia Reference Référence	Capacidad de producción (mín./máx.) Production capacity (min./max.) Capacité de production (min./max.) l	Dimensiones útiles cámara Usable chamber dimensions Dimensions utile chambre Ø x H mm	EXTERIOR L x D x H mm	Temp. °C	Presión máx. Max. pressure Pression max. Barg	Potencia ¹ Power Puissance W	Peso bruto Gross weight Poids brut Kg	Tensión ² Voltage Tension V	Frecuencia Frequency Fréquence Hz
AE-20-MP³	2/18	320 x 350	615 x 815 x 735	50-125	1,4	3000	95	230	50/60
AE-40-MP	3/37	320 x 500	615 x 815 x 1100	50-125	1,4	6000	165	400	50/60
AE-60-MP	3/57	320 x 700	615 x 815 x 1320	50-125	1,4	9000	180	400	50/60
AE-80-MP	5/78	420 x 500	755 x 935 x 1285	50-125	1,4	15000	250	400	50/60
AE-100-MP	5/97	420 x 700	755 x 935 x 1385	50-125	1,4	15000	295	400	50/60

1. Modelos especiales con potencia incrementada disponibles bajo pedido, consultar opciones / Special models with increased power available under request, consult for options / Modèles spéciaux avec puissance accrue disponibles sur demande, consulter les options

2. Consultar por otras tensiones / Ask for other voltages / Demander d'autres tensions

3. Incluye mesa de acero inoxidable con ruedas (+info: en Accesorios) / Included stainless steel table with casters, (+info: in Accessories) / Table en acier inoxydable incluse avec roulettes (+info: dans Accessoires)

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Intended use

• Automation and acceleration of the dispensing phase in repetitive operations of medium and small volumes controlled by optical sensor or pedal control.

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Components supplied

- 1 Control box with timer.
- 1 Dosing foot pedal.
- 1 Support with proximity sensor (for object detection, for example: vessel).
- 2 Clamps for tube.
- 1 Dosing tube with nozzle.
- 1 Power cord.
- 1 Sensor connection cord.
- 1 Manual of instructions.

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Referencia: **PH-MP**

PH METER

- Control and registry of pH, with real-time monitoring on screen.

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Referencia: **KLL-MP**

AUTOMATIC WATER FILLING KIT

- Water pump to automate the feed of the water tank with purified water in media preparators AE-MP series.
- Compatible with installations that either have a purified water network, a purified water tank or facilities that have an unpurified water network, in the latter case, the kit should be supplied with two other accessories: water purifier (ECOPUR-500) and purified water tank (TANK-KLL).
- Must be installed in our facilities.

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Referencia: **CAP-MP**

EXTERNAL PROBE ADAPTER

- Plug with special adapter for an external rigid probe for continuous validation processes.

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Referencia: **SW8000**

PC CONNECTION SW8000

- Communication software between the equipment and the PC.
- Used with Ethernet connection.
- USB adapter.
- Compatible with touchscreen circuits.

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Referencia: **TABLE-MP**

TABLE

- Stainless steel table with wheels (with brakes on two of them).
- This table is included with the AE-20-MP model.
- Dimensions (LxDxH): 700 x 700 x 600 mm.

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Referencia: **CP-MP**

COMPRESSED AIR SYSTEM

- Air compressor that offers support pressure during cooling phase to avoid liquid loss by evaporation. This accessory is necessary to enable external dispensation control through the Dosing station for culture media (DW-MP).
- Must be installed in our facilities.

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Intended use

- Incineration of organic matter to prevent cross contamination by contact or splashing.
- Specially indicated in microbiology for the sterilization of small metal utensils like needles, lancets, inoculating loops and tweezers.

Features

- Quartz incinerator tube, with electric element and thermally insulated with ceramic fiber.
- Protective external grid made of stainless steel. Epoxy resin oven painted base.
- Start up pilot light.
- Sterilization by infrared rays.
- Organic material is incinerated by introducing it into an incandescent quartz tube, avoiding splashing of infectious microorganisms.

Advantages

- Perfect substitute of Bunsen burners and alcohol lamps.
- Sterilization is completed within 5-7 seconds at 850 °C.
- Very safe to operate with, no open flame while functioning.

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Kjeldahl

The Kjeldahl method is a chemical analysis process to determine nitrogen content. It is widely used as a reference method to estimate protein content in food.

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Randall/Twisselman

The Randall/Twisselman methods are used to extract substances from a solid sample, usually of lipidic nature, through a solvent.

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Weende

The Weende method is commonly used to quantify crude fibre in a food sample after an acidic and alkaline hydrolysis treatment.

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Van Soest

The Van Soest method is used to determine acid detergent fibre (ADF) and neutral detergent fibre (NDF) content in a sample based on the indigestible residue obtained.

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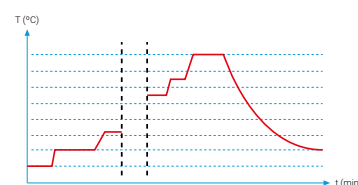
Intended use

- Analysis of proteic nitrogen in organic substances such as nourishing products, fertilizers, seeds, plants, beer, meat, sauces, water or petroleum. See page 66.
- Sample pretreatment (hydrolysis) for fat extraction.
- Soil digestion for metal analysis.

Advantages

- Programmable autostart.
- Built-in temperature regulator which has 9 programs with 18 segments of temperature, ramp and time each.

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- Test tube rack with support (necessary tubes included).
- Fume collector with support.

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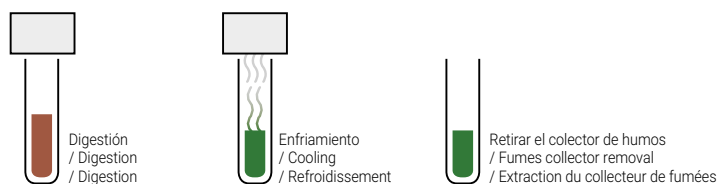
Features

- Quick and safe manipulation of samples.
- Integrated support for sample tubes enabling an easier cooling and saves space.
- Upper chassis made of aluminum alloy and Halar® coating. Metallic bottom chassis painted with epoxy resin. Metal block made of aluminum alloy.
- Thermal insulation: rock wool and ceramic fiber.
- Safety thermostat to prevent overheating.
- Contains calibration port for external probes.
- Ideal for use in combination with the **SCRUBBER Fumes neutralization unit**, (see page 70). Both systems are connected by a flexible tube that joins the Compact digestion system fume collector (MBC) and the Fumes neutralization unit (Scrubber).

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Intended use

- Extraction and neutralization of harmful fumes.

Features

- The gas produced during the digestion is extracted by the action of a vacuum pump. First, the gas undergoes a phase of condensation that acts as preliminary extractor for steams and dragged liquids, avoiding warming or volume increase in posterior wash-up solution.
- The acid or alkaline vapors are washed and neutralized in the next step. In this neutralization phase, most of the unwanted particles are retained through granules of activated charcoal. The used air in the process is conducted to an extractor or to the exterior.
- Case made of stainless steel AISI 304 painted with epoxy resin. Glass objects made of: Borosilicate 3.3. Pipes made of: PVC (Cristalflex®), silicone and Teflon®. Pump made of: PPS, EPDM and Viton®.
- Ideal to use together with **MBC Compact digestion system** (see page 68). Both systems are connected by a flexible tube that joins the Compact digestion system fume collector (MBC) and the Fumes neutralization unit (Scrubber).

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Advantages

- Vacuum pump with acoustic insulation of 48l/min.
- Filtration and neutralization of harmful fumes by a filter of activated carbon.
- Absorption valve manually adjustable from 10 mbar to 800 mbar.
- Adjustable steam suction speed.
- Closed fumes evacuation system.

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Intended use

• Distillations for the Kjeldahl analysis, ammoniacal nitrogen, alcoholic strength by volume, phenols, volatile acids and cyanides.

Features

- System control by microprocessor with LCD screen.
- Additionally, the system has fixed programs:
 - Ammonium sulphate test.
 - Preheating of the system.
 - Sample waste washing.
- Adaptable system for distillation tubes of different diameters.

Advantages

- Kit for automatic titration included for DNP-3000 and optional for DNP-2000MP.
- Programs: 11 predefined (editable) and 9 free.
- Calibrable pumps by the user.
- Steam generator with water level control.
- Sample suction (only for DNP-2000MP/3000).
- Boric acid pump (only for DNP-2000MP/3000).
- Results in N or total protein (only for DNP-3000).

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Safety

- Open door sensor.
- Distillation tube detection.
- Over-temperature thermostat.
- Overpressure pressure switch.
- Pressure regulator at the cooling water inlet.

Performance

- Measurement range: from 0.1 to 200 mg of N.
- Nitrogen recovery: $\geq 99.5\%$.
- Reproducibility: $\pm 1\%$.
- Distillation time: 4 min/100 ml.
- Electric power: 2000 W (100%).

Programmable parameters and values

- Water for dilution: 0-240 ml.
- NaOH solution: 0-240 ml.
- H_3BO_3 solution: 0-240 ml (only for DNP-2000MP/3000).
- Delay time: 0-30 min.
- Distillation time: 0-30 min.
- Steam flow regulation: 30-100%.

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- 1 distillation tube Ø42 x 300 mm.
- 2 fast connection H_2O and NaOH containers of 10 litres.
- 1 fast connection H_3BO_3 container of 5 litres (only for DNP-2000MP/3000).
- Additional pipe without filter for steam addition (only for DNP-2000MP/3000).
- Automatic titrator (only for DNP-3000).
- Several connection hoses.
- Anti-drip tray.

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Intended use

- Quantification of nitrogen and protein.

It includes

- Magnetic stirring, reaction vessel and complete titrator.

Features

- Potentiometric titrator.
- Multiple titration programs.
- pH calibration.
- Burette autocalibration.
- Results in N or total protein.
- 2 USB ports:
 - 1 to copy tests and extract data.
 - 1 to connect printer, keyboard or mouse.
- Ethernet port: optional software connection.

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Referencia: **SOFT-TITRA**

PC CONNECTION SOFT-TITRA

- Communication software between the titrator and the PC.
- Connection via Ethernet.
- Control to start and stop analysis.
- Show real-time data directly from the workstation.
- Manage data stored locally or on a server (search, compare, delete, print, etc.).
- Export data.

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Referencia: **ADAP-BU**

DNP ADAPTER FOR BÜCHI® TUBES

- Adapter to use the Büchi® tubes in the DNP distiller.

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Intended use

• Extraction system for fat or other soluble materials according to Randall or Twisselman methods, based on the evolution of the traditional Soxhlet method. For the analysis of food, feed, soil, detergents, polymers, paper pulp, fibers, gums, textiles, pharmaceuticals and petrochemicals.

Features

- Heating by an electric plate with armoured heater to assure homogeneous distribution of heat.
- Control of temperature by P.I.D. microprocessor action and probe Pt100 class A.
- Over temperature safety control by independent thermocouple.
- Vitón® sealing.
- In accordance with standardized international methods: AOAC, ISO, AACC.
- Case made of stainless steel AISI 304 painted with epoxy resin.

Advantages

- Independent programming of temperature and extraction times for "boiling", "rinsing" and "recovery" phases.

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Technical Data

- Temperature range: room temp. +5 °C to 220 °C.
- Solvent volume: up to 50 ml.
- Solvent recovery: up to 95%.
- Reproducibility: ±1%.
- Extraction thimbles: Øext 26 x 60 mm.
- Programmable boiling extraction time.
- Programmable rinsing extraction time.
- Programmable solvent recovery time.
- Grade of protection: IP65.

Components supplied with the equipment

- 12 extraction vessels + magnetic adapters.
- 6 thimble supports.
- 2 racks for extraction cartridges.
- 2 racks for extraction vessels.
- Aligning rack.
- Insertion handle.
- 2 silicone tubes 1,5m (Ø8x14).
- Cellulose thimbles box (25 units).
- Vessel manipulation rack.
- 6 EPDM seal set.

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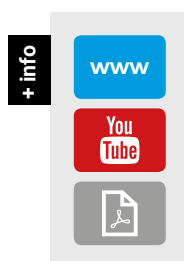
Intended use

• Determination of raw/crude fibre content by *Weende*, acid detergent fibre (ADF) and neutral detergent fibre (NDF) by *Van Soest* and acid detergent lignin (ADL).

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Features

- Extraction and filtration without sample transfer.
- High reproducibility of results.
- Up to 36 analysis/day (*Weende* method).
- Versatile use: samples can be dried/weighted at every phase if needed.
- System for 6 simultaneous samples.
- All electric and mechanic elements are conveniently protected.
- Case made of stainless steel AISI 304 painted with epoxy resin.
- Additionally, it can be used in combination with the **Cold fat extraction system EF-6P** (see page 82), to guarantee the optimum working conditions of the system.

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Technical Data

- Range of measurement: 0.1-100%.
- Reproducibility: $\pm 1\%$ for samples with a fiber content ranging from 1-30%.

Components supplied with the equipment

- 6 glass crucibles porosity P-2.
- Crucible rack.
- Heater lid.
- Crucible handling support.
- Crucible vessel tong.
- Connection hoses.

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Pressure contribution

- F-6P has a peristaltic pump that provides the necessary conditions for filtering.

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Intended use

- Designed for cold fat extraction with acetone or other solvents.
- Preliminary degrease before hot fibre extraction of samples with more than 1% fat content.

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Features

- Up to 6 simultaneous samples per cycle.
- Case made of stainless steel AISI 304 painted with epoxy resin.
- Ideal for use in combination with the **Manual fibre determination system F-6P** (see page 80).

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Pressure contribution

- EF-6P has a peristaltic pump that provides the necessary conditions for filtering.

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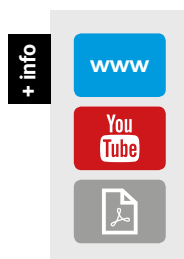
Intended use

• Determination of the alcoholic strength by volume in wines, musts, beers and spirit drinks of high graduation according to the regulation OIV, CEE 2676/90 and CE 2870/2000. Volatile Acidity and Sorbic Acid can also be measured.

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Features

- Distillation by steam dragging.
- Steam generator automatically fed with distilled water, having electronic level control and automatic stop in the case of lack of water in the container.
- Safety thermostat.
- Automatic distillation stop.
- Manual levelling of the distillate avoiding adding distilled water.
- Automatic stop of refrigerating water at the end of distillation.
- Optical display of minimal level of water in the steam generator.
- Button to add distillate manually.
- Case made of stainless steel AISI 304 painted with epoxy resin.

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Technical data

- Distilled water reservoir: 10 l.
- Approx. distillation time: 7 min./200 ml.

Components supplied with the equipment

- 1 distillation tube for maximum sample volume of 100 ml.
- 1 distillation tube for maximum sample volume of 250 ml.
- 1 volumetric flask of 100 ml.
- 1 volumetric flask of 200 ml.
- 1 volumetric flask of 250 ml.
- 1 fast connection container of 10 litres for H₂O.
- Connection hoses.

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TB-42300 TB-80300

GRADILLA 6 TUBOS / 6 TUBES RACK / PORTOIR 6 TUBES

Referencia Reference Référence	Ø x H mm
GRA-640	Ø 42 x 300
GRA-680	Ø 80 x 300

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Technical features

- Intuitive alphanumeric user interface through a 3,5" LCD touchscreen that allows the display of time, temperature and cycle evolution in graphic format in real time.
- Compatible with Celsius or Fahrenheit scales with a resolution of 0.1°C/°F.
- Capable of advanced cycle and timer programming, with up to 10 preset programs editable in parameters and name, programmable at a fixed temperature during an infinite time or with up to 10 time and temperature segments. Timer start and timer stop customizable by date and time.
- Audiovisual security alarms for cycle finalization, probe failure, heating error, overheating and temperature out of range.

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Technical features

- All incubation and drying equipment have a Ethernet connection that allows the use of the RAYPANet online platform on a PC to view and export all the information of the cycles with graphs and all the relevant data, with the ability to export data reports.
- All cycle data can be viewed on the platform in real time or through the cycle history registry. Furthermore, all data can be exported in .csv and .pdf file formats for later consultation, study and storage.
- The platform is compatible with all web browsers and multiple equipment can be connected and controlled simultaneously.

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Intended use

• Incubation of microbiological and cell cultures, crystal growth, conditioning of thermosensitive materials and controlled storage of samples.

Features

- Incubation by natural convection.
- Temperature range from room temperature +5°C to 80°C.
- Temperature and time controlled by digital PID microprocessor, with 10 editable programs of customizable name, programmable with up to 10 temperature/time segments.
- Intuitive alphanumeric touchscreen user interface.
- Pt100 class A temperature probe.
- Ethernet connection.
- External metal frame with epoxy resin coating.
- Inner frame, dual chamber, inner side of external door and perforated trays made of stainless steel AISI-304.
- Tempered glass inner door with magnetic lock.
- Silicone rubber door seal.
- Only for ID-480 and ID-720: Double-leaf inner and outer doors with magnetic lock and 4 casters with brakes.

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Safety

- Audiovisual security alarms.
- Safety thermostat.
- Stainless steel interior and corrosion-resistant, easy to clean metal exterior.
- Homogeneity and temperature stability according to DIN 12880 regulation.
- Safety according to EN 61010-1 and EN 61010-2-010 regulations.

Supplied with

- 2 perforated trays made of stainless steel AISI-304 adjustable in height, with anti-tilting system.
- Access to the online platform RAYPanet.

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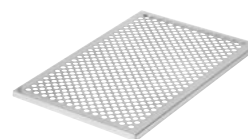
Referencia: **FIL-ID/DOD**

BACTERIOLOGICAL FILTER

- Bacteriological filter of 0.22µm for inlet and/or exhaust air.

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Referencia: **MCLL**

LOCKABLE HANDLE

- Lockable door protects samples from unauthorized access.

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GLASS DOOR

- Double tempered glass exterior door in special models.

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Intended use

• Water analysis, B.O.D. test (Biological Oxygen Demand), enzymatic assays, cell and bacteria culture at low temperature, botanics, aging, cosmetics and sample storage.

Features

- Forced air cooling incubation.
- Temperature range from 0°C to 45°C.
- Temperature and time controlled by digital PID microprocessor, with 10 editable programs of customizable name, programmable with up to 10 temperature/time segments.
- Intuitive alphanumeric touchscreen user interface.
- Pt100 class A temperature probe.
- Ethernet connection.
- Two free plugs inside.
- Internal illumination.
- External probe access.

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Safety

- Audiovisual security alarms.
- Safety thermostat.
- Corrosion resistant and easy to clean powder coated metallic exterior
- Safety according to EN 61010-1 and EN 61010-2-010 regulations.

Supplied with

- 3 removable plastic-coated stainless steel trays (4 trays for model IRE-475).
- Access to the online platform RAYPAnet.

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Intended use

• Drying and sterilization of samples in all types of research and industrial environments.

Features

- Drying by natural convection.
- Temperature range from room temperature +5°C to 250°C.
- Temperature and time controlled by digital PID microprocessor, with 10 editable programs of customizable name, programmable with up to 10 temperature/time segments.
- Intuitive alphanumeric touchscreen user interface.
- Pt100 class A temperature probe.
- Ethernet connection.
- External metal frame with epoxy resin coating.
- Inner frame, dual chamber, inner side of door and perforated trays made of stainless steel AISI-304.
- Silicone rubber door seal.

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Supplied with

- Two perforated trays made of stainless steel AISI-304 adjustable in height, with anti-tilting system.
- Access to the online platform RAYPanet.

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Intended use

• Drying and sterilization of samples in industrial and research environments in the chemical and technological sector such as drying of varnishes and solvent coatings that require standardized, safe and reproducible drying processes.

Features

- Forced air drying.
- Temperature range from room temperature +5°C to 250°C.
- Temperature and time controlled by digital PID microprocessor, with 10 editable programs of customizable name, programmable with up to 10 temperature/time segments.
- Intuitive alphanumeric touchscreen user interface.
- Pt100 class A temperature probe.
- Ethernet connection.
- External metal frame with epoxy resin coating.
- Inner frame, dual chamber and perforated trays made of stainless steel AISI-304.
- Double insulated walls and door with low conductivity glass wool.
- Silicone rubber door seal.
- Only for DAF-425 and DAF-635: Double-leaf door and 4 casters with brakes.

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Supplied with

- 2 perforated trays made of stainless steel AISI-304 adjustable in height, with anti-tilting system.
- Access to the online platform RAYPANet.

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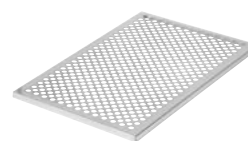
Referencia: **FIL-DAF**

BACTERIOLOGICAL FILTER

- Bacteriological filter of 0.22µm for inlet and exhaust air.

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Intended use

• Materials that require effective drying and are sensitive to heat such as electronic components, food, small components, pharmaceuticals and chemicals.

Features

- Vacuum drying.
- Temperature range from +35°C to 250°C.
- Temperature and time controlled by digital PID microprocessor, with 10 editable programs of customizable name, programmable with up to 10 temperature/time segments.
- Intuitive alphanumeric touchscreen user interface.
- Pt100 class A temperature probe.
- Ethernet connection.
- Inert gas connection.
- External metal frame with epoxy resin coating.
- Inner frame and dual chamber made of stainless steel AISI-304.
- Reinforced double glass door.
- Double insulated walls and door with low conductivity glass wool.
- Galvanized aluminum perforated trays.
- Silicone rubber door seal.

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Safety

- Safety double glass against implosions.
- Audiovisual security alarms: end of cycle, probe failure, heating error, overheating and temperature out of range.
- Safety thermostat.
- Stainless steel interior and corrosion-resistant, easy to clean metal exterior that remains at safe temperature without overheating.
- Safety according to EN 61010-1 and EN 61010-2-010 regulations.

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Supplied with

- 2 perforated removable trays made of aluminium with galvanic treatment.
- Access to the online platform RAYPANet.

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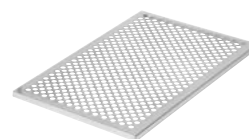
LOCKABLE HANDLE

- Lockable door protects samples from unauthorized access.
- Reference: **MCLL**

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VACUUM GENERATION SET FOR EV-50

- System of 3 parts: settling vessel, filter and vacuum pump.
- Includes 2 meters of silicone tube (Ø 8 x 14 mm).

Reference: **VGS**

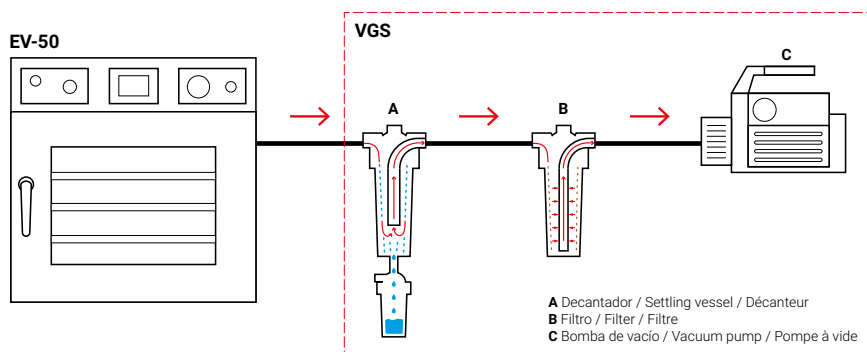
- **Clean room** VGS adaptation available under special request.

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- Settling vessel with large storage capacity.
- Fastening system by central rod for attachment and reliable sealing.
- Emptying system without interruption of the vacuum supply (settling vessel).

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- Filtering elements with a filtration capacity of 0.1 micron.
- Fastening system by central rod for attachment and reliable sealing.

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- Two-stage vacuum pump with vacuum gauge and non-return valve to protect the system from oil contamination.
- Ergonomic handle with protection, for a more comfortable grip.
- Visor with oil level indicator and oil drain system for easy emptying.

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Intended use

- Incubation of water samples, reagent heating, substrate fusion, cell culture incubation, etc.
- Compatible with water or oil.

Features

- Compatible with water and oil with a temperature range of + 5°C ambient temperature to 200°C.
- Temperature controlled by digital PID microprocessor, with 10 editable programs of customizable name, programmable with up to 10 temperature/time segments or at a fixed temperature.
- Timer start and timer stop adjustable by minute, hour, day, month and year.
- Intuitive alphanumeric touchscreen user interface.
- Pt100 class A temperature probe.
- Inner tank and external case made of AISI-304 stainless steel.
- Heating element inside the tank with a protecting grid.
- Incorporated drainage tap.
- Maximum and minimum tank filling: 50% - 80% of total capacity.

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Advantages

- Possibility to select between Celsius or Fahrenheit degrees.
- 3.5" color touch screen.
- Graph visualization: programmed objectives and real curve.
- Visualization of program phases.

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Safety

- Visual and acoustic alarms: probe failure, heating failure, temperature outside the chosen range (both over and under).
- Safety under regulation EN 61010-1.
- Safety thermostat with manual rearms according to DIN 12877 class 2.

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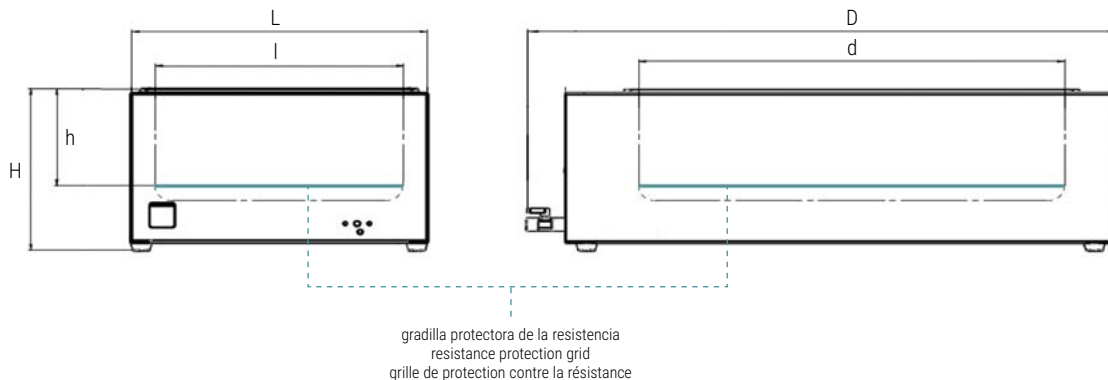
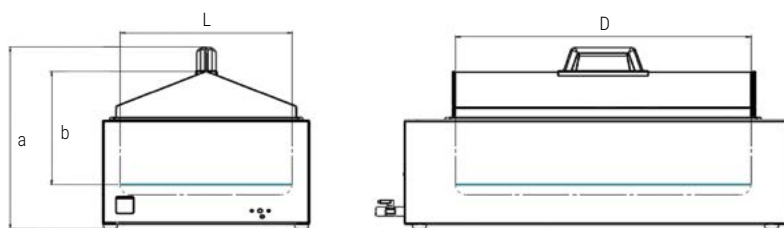
MARCO SOPORTE * / SUPPORT FRAME / SUPPORT PORTOIR

Referencia Reference Référence	Serie baño Bath series Série bain
MB001	WBD-5
MB002	WBD-12
MB003	WBD-20
MB003 (x2)	WBD-40

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Intended use

- For cases where homogeneity is required at high temperatures with gentle heating.
- Indicated for ovoid materials in drying operations, evacuation, etc.

Features

- Adjustable temperatures on plate surface up to 400 °C.
- Temperature regulation by energy impulses regulator.
- Steel hot plate with surrounding heating elements.
- Stainless steel upper case.
- Control panel with pilot-lights indicating connection to the mains and heater functioning.
- Removable stainless steel inner tank, enabling to use of the apparatus as a heating plate.

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Follow us:

ONLY RAYPA MEDICAL AUTOCLAVES COMPLY WITH ISO 13485:2016 CERTIFICATION

RAYPA

Leading
Lab Technologies

Avinguda del Vallès, 322
Pol. Ind. "Els Bellots"
08227 Terrassa (Barcelona) Spain

raypa@raypa.com
www.raypa.com

Tel. +34 937 830 720

R. ESPINAR, S.L.